

# SIMPLIFIED ROUTINE QC MONITORING OF GLYCANS USING THE ACQUITY™ QDA™ II MASS DETECTOR

Henry Foddy<sup>1</sup>, **Scott Berger**<sup>2</sup>, David Douce<sup>1</sup>, Cristian Cojocariu<sup>1</sup>, Ali Al Kaabi<sup>3</sup>, Andreas Fluetsch<sup>3</sup>

<sup>1</sup>) Waters Corporation, Wilmslow, United Kingdom

<sup>2</sup>) Waters Corporation, Milford, MA

<sup>3</sup>) Johnson & Johnson Innovative Medicine, Bern, Switzerland

## INTRODUCTION

Glycans are complex carbohydrate structures that, when covalently linked to proteins and lipids, play an integral role in therapeutic and diagnostic processes. Precise and reliable glycan profile analysis is essential, often requiring comprehensive approaches to elucidate both structural and functional properties.

Typically, these workflows involve multi-step processes that leverage advanced mass spectrometry techniques, commonly used for their selectivity and sensitivity advantages. However, such techniques are often associated with high operational costs and complexity, which can make routine monitoring challenging. Consequently, there is a growing interest in exploring cost-effective alternatives for these workflows, such as the ACQUITY QDa II Mass Detector, that may be better suited to high throughput, robust applications, particularly in QC testing laboratories.



Figure 1. The ACQUITY QDa II Mass Detector and ACQUITY Premier UPLC™ configuration

A cost-effective and reliable instrument, the ACQUITY QDa II Mass Detector offers an extended mass range of 30–1500  $m/z$ , as well as the sensitivity and robustness required for such routine pharmaceutical applications. Automated resolution, calibration, and tuning on start-up allow for simplicity and ease-of-use, reducing the burden on the analyst and facilitating more sustained instrument uptime.

**TO DOWNLOAD A COPY OF THIS POSTER, VISIT [WWW.WATERS.COM/POSTERS](http://WWW.WATERS.COM/POSTERS)**

In this study, a glycan fragment was monitored and quantified using the ACQUITY QDa II Mass Detector, using intentional in-source fragmentation to target a fragment ion with a mass of 844 Da. This targeted fragmentation approach offers a unique and reliable method for fingerprinting and monitoring larger glycan structures, making it well-suited for routine QC testing in a cost-effective manner. The workflow was evaluated for sensitivity, linearity, compound carryover, and overall method performance.

## RESULTS & DISCUSSION

A method was developed using an unnamed 20 µg/mL glycan sample. The instrument was operated in positive electrospray ionization (ESI+) mode with an optimized cone voltage of 35 V, a capillary voltage of 1.2 kV, and a desolvation temperature of 600 °C. The fragment ion of interest was monitored at  $m/z$  845 using selected ion recording (SIR) mode. Simultaneously, a full scan acquisition (total ion current, TIC) from  $m/z$  150 to 1500 was performed to evaluate the overall elution profile of the protein structure. An example of the data from these two monitoring channels is shown in Figure 2

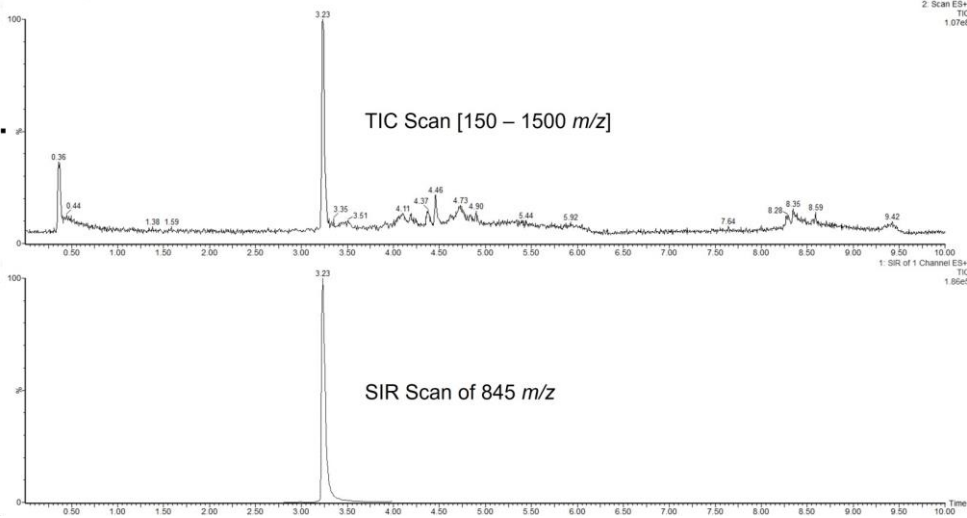


Figure 2. Stacked chromatogram plots showing a TIC Scan from  $m/z$  150 – 1500 and an SIR channel monitoring  $m/z$  845 in positive (+ve) electrospray ionization mode, for the same injection of a 20 µg/mL glycan sample

### Carry-over Assessment

An ACQUITY Premier Protein BEH™ C4 300 Å 1.7 µm, 2.1 x 50 mm column was selected for the work, and a 30 °C column temperature was found to be optimal using a 1 µL injection. An ACQUITY Premier LC System was also used in this workflow, with the FTN (flow-through needle) design helping to reduce sample dispersion and carryover. Reducing carry-over was a critical aspect of this work. To address this, fragment ion carry-over was assessed, with the method modified to minimize carry-over in subsequent blank injections.

Incorporating Isopropyl alcohol (IPA) into the composition of the mobile phase was pivotal in this solution, reducing carry-over in solvent injections immediately following the sample to less than 1%, while maintaining the instrument's response to the compound (see Figure 3).

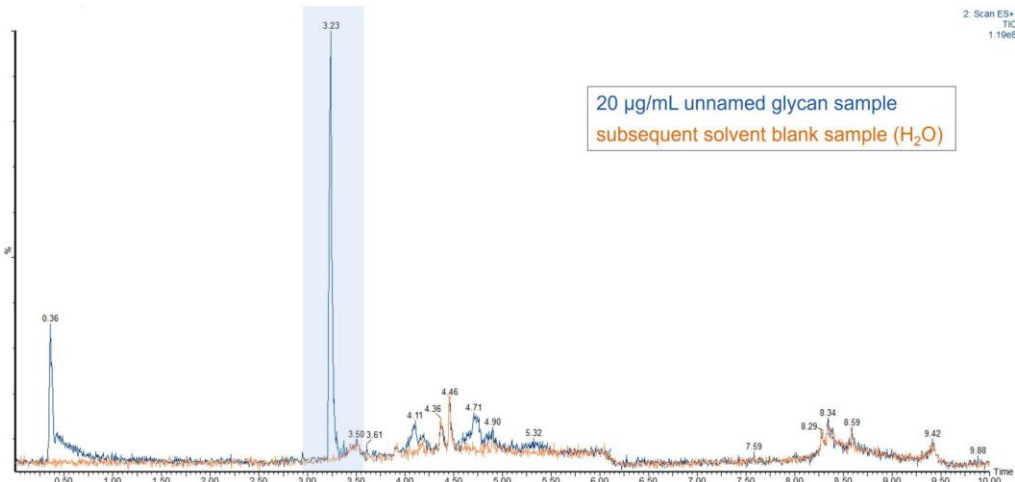


Figure 3. Overlaid TIC chromatograms of a 20 µg/mL injection of an unnamed glycan sample (blue) and the subsequent solvent blank injection (orange)

## RESULTS & DISCUSSION

### Linearity

The peak area response of the fragment ion of  $m/z$  845 was linear over a calibration range of 2 - 20 µg/mL (corresponding to 2 - 20 ng of glycan sample on column) with a coefficient of determination  $R^2 = 0.99$ . No internal standard correction was used. The corresponding residual values for any given datapoint did not exceed 12%, as shown in Figure 4. Precise and reliable quantitation offers the possibility of accurate routine determination of this fragment ion.

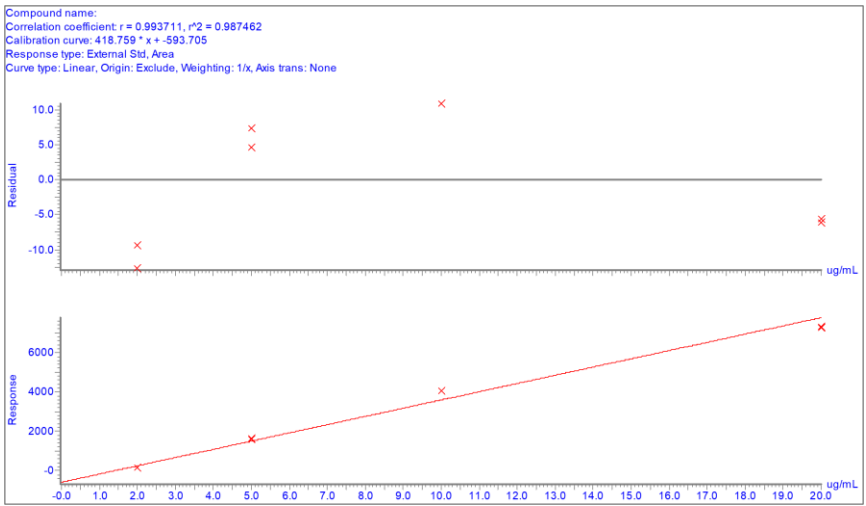


Figure 4. Calibration response of the  $m/z$  845 fragment ion over a concentration range of 2 – 20 µg/mL. Residual values (top) as well as calibration curve (bottom) are shown.

### Quantitation

A sample was also assessed at 1 µg/mL (corresponding to 1 ng on-column) giving a signal-to-noise ratio of 102:1 (S/N as peak to peak) which can be seen in Figure 5.

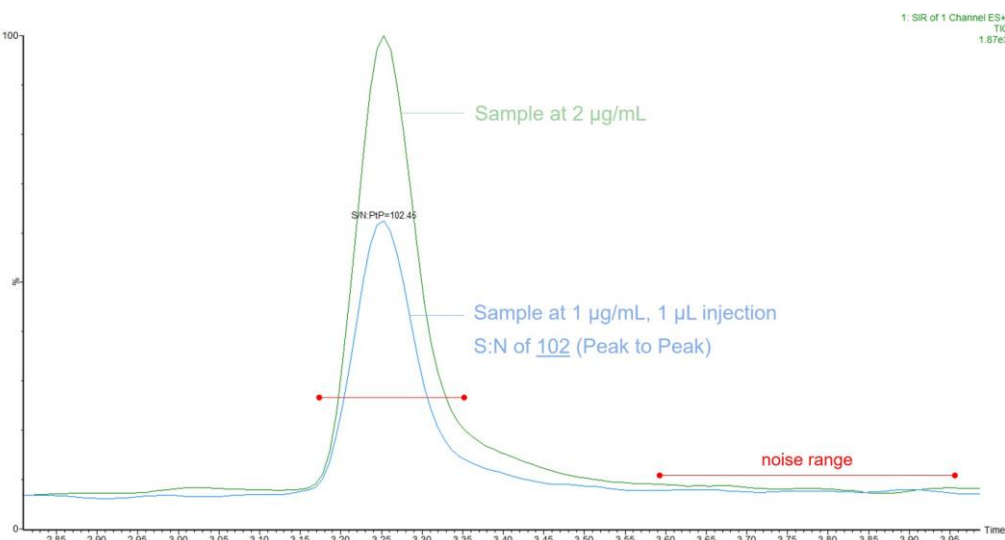


Figure 5. An SIR ( $m/z$  845) chromatogram of a 1 µg/mL glycan sample with a S/N (peak to peak) of 102, with signal and noise regions selected shown in red

## CONCLUSION

The ACQUITY QDa II Mass Detector is a compact and stackable device designed for seamless integration with LC systems. It is user-friendly, robust, and cost-effective, making it an ideal choice for routine glycan monitoring. It offers:

- Adequate sensitivity, detecting glycan fragments ions at 1 ng on-column in challenging matrices
- Robust performance, handling challenging matrices with reduced carryover and ensuring accurate quantitation
- Linear response for glycan fragment ion over a concentration range of 2 to 20 µg/mL
- High-throughput capability to support efficient quality control for complex proteins, reducing the need for more expensive equipment
- Reduced carryover and accurate quantitation of challenging matrices by employing the ACQUITY Premier Protein BEH C4 Column, and ACQUITY Premier LC System
- Compatibility with both MassLynx™ and Empower™ Software for seamless integration into QC labs

ACQUITY, QDa, UPLC, BEH, MassLynx, and Empower are all trademarks of Waters Corporation