

Automated Cyclic Ion Mobility–Mass Spectrometry MS/MS Workflow and Data Analysis for Shotgun Lipidomics

Ballal Hossen¹, Nyasha Munjoma², Susan Slade², Emma Marsden-Edwards², Johannes PC Vissers^{2*}, Dominik Schwudke^{1,3,4*}

¹ Division of Bioanalytical Chemistry, Research Center Borstel - Leibniz Lung Center, Borstel, Germany, ² Waters Corporation, Wilmslow, SK9 4AX, United Kingdom, ³ German Center for Infection Research, Thematic Translational Unit Tuberculosis, Borstel, Germany, ⁴ German Center for Lung Research (DZL), Airway Research Center North (ARCN), Borstel, Germany

Introduction

Recent advances in technology enable sensitive lipid identification down to the single-cell level. Improved MS techniques provide greater precision in characterizing cell-cell interactions, cellular states, and phenotypes based on lipidomic profiles. However, even state-of-the-art lipidomics approaches remain limited in their ability to address the heterogeneity and complexity of cellular lipidomes. Here, the integration of cyclic IM into a shotgun lipidomics workflow is described. In proof-of-principle experiments, this additional dimension of separation improves spectral clarity and facilitates the resolution of isobars and isomers beyond m/z -based discrimination. An automated acquisition and data processing workflow was developed to handle two-dimensional data (m/z and drift time) for both MS1 and MS2 of quadrupole-selected precursors.

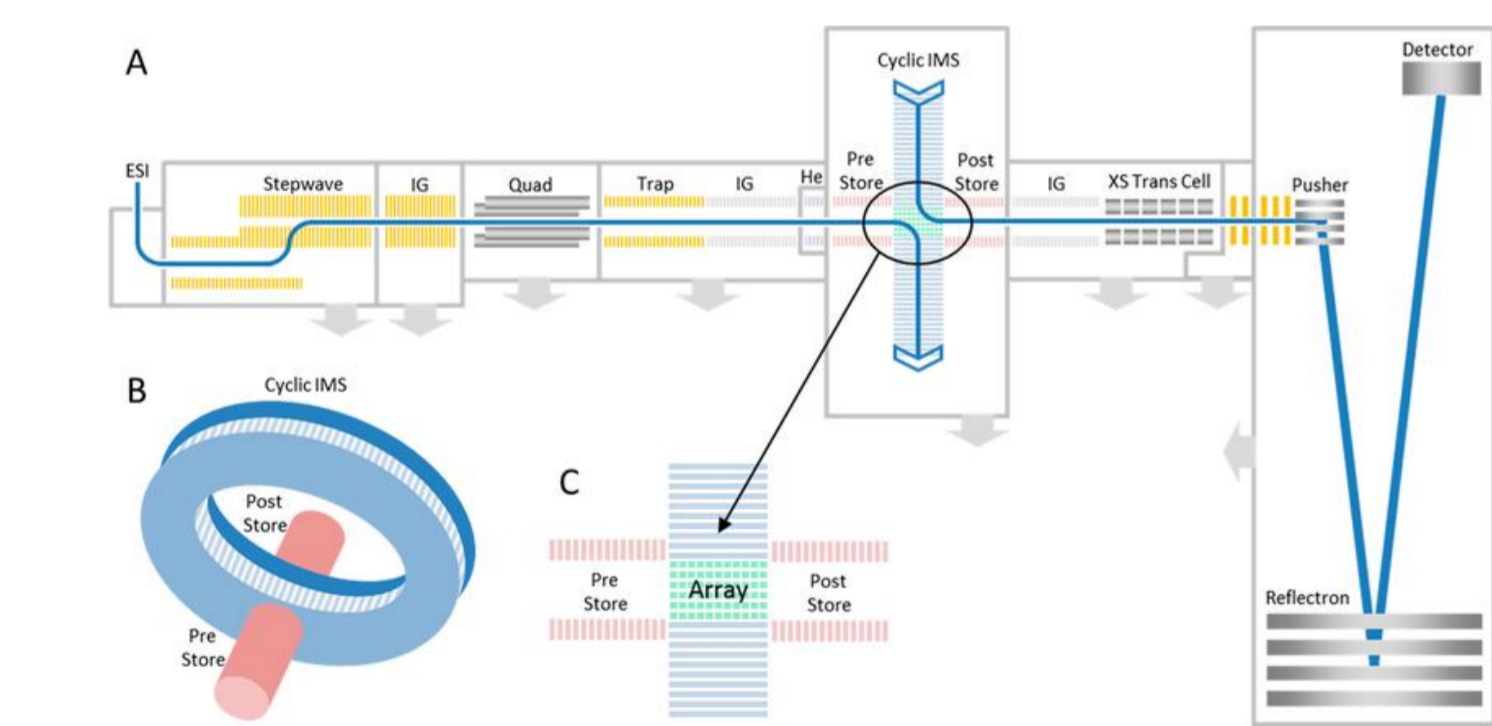


Figure 1. SELECT SERIES™ Cyclic™ IMS [1]. (A) Overall schematic, (B) and (C) cIM array.

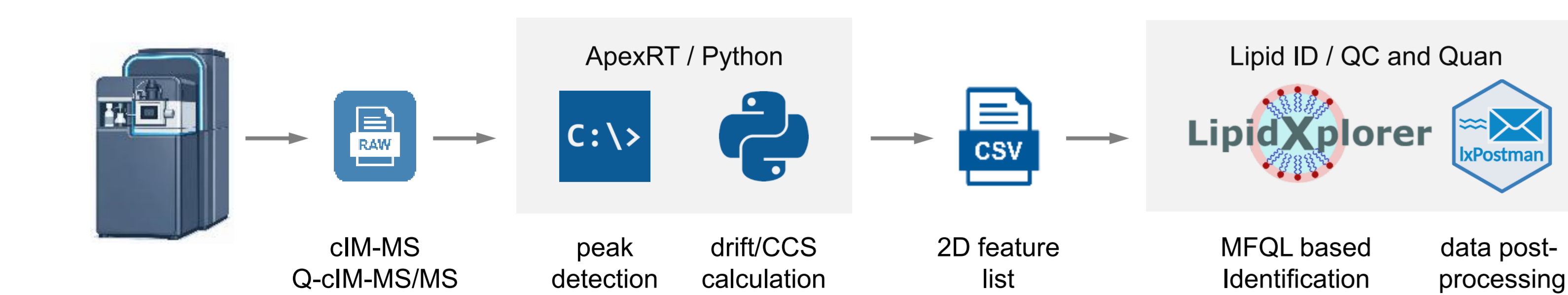


Figure 2. Semi-supervised prototype data acquisition and processing workflow, illustrating multidimensional peak detection and lipid annotation/identification, using development software tools, LipidXplorer [3] and IxPostman [4].

Experimental

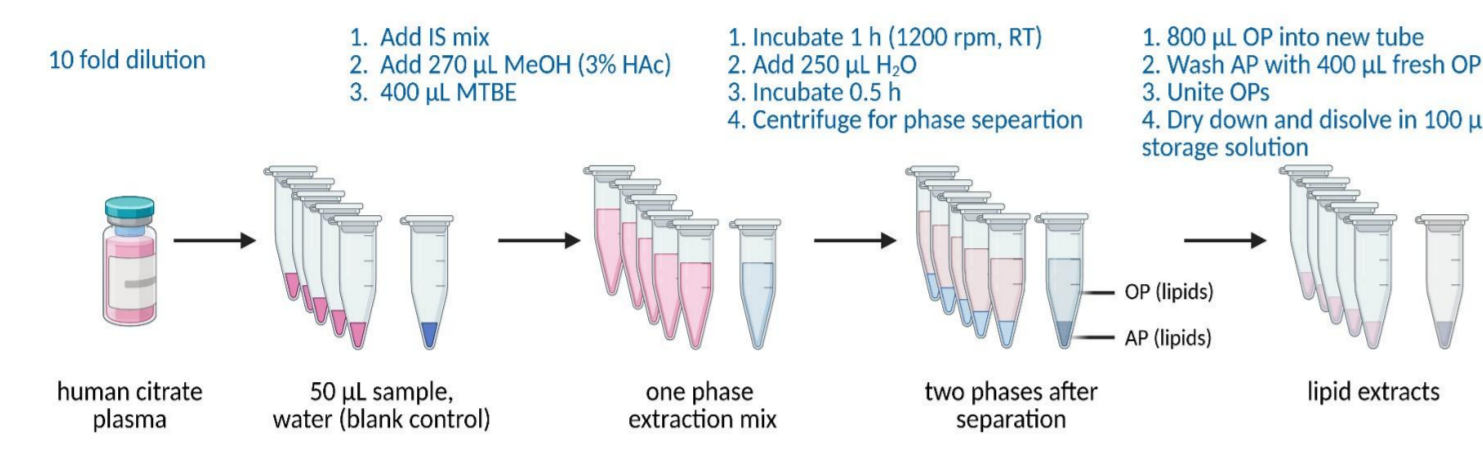


Figure 3. Sample preparation for shotgun lipidomics. Five independent lipid extractions were performed using a customized MTBE protocol [2]. Extracts are dissolved in MS-mix in 1/20 dilution directly before measurements.

Created in BioRender. Schwudke, D. (2026) <https://BioRender.com/fswd2xa>

Direct Infusion Mass Spectrometry

- SELECT SERIES Cyclic IMS (Waters Corporation)
- nano-ESI (+) using Triversa NanoMate (Advion) 1.1 kV at 1.0 Psi
- Source temp: 150 °C
- Acquisition methods: HDMS (cIM-MS), HDMS/MS (Q-cIM-MS/MS) in v-mode

Data processing

- MassLynx™ v4.2 and DriftScope™ 3.0 (data acquisition/review)
- ApexRT 1.17 (peak detection)
- Pandas, numpy Python libraries (drift time calculation)
- Prototype software LipidXplorer 1.5 [3] and IxPostman [4]

Results

Experiment were designed using Cyclic IMS Method Editor, Figure 4. A representative direct-infusion cIM-MS 2D map is shown in Figures 5A and 5B. These data demonstrate the ability of cIM to resolve isobaric, Q-isolated precursor ions, resulting in reduced spectral complexity. Results are shown in Figures 6 to 9.

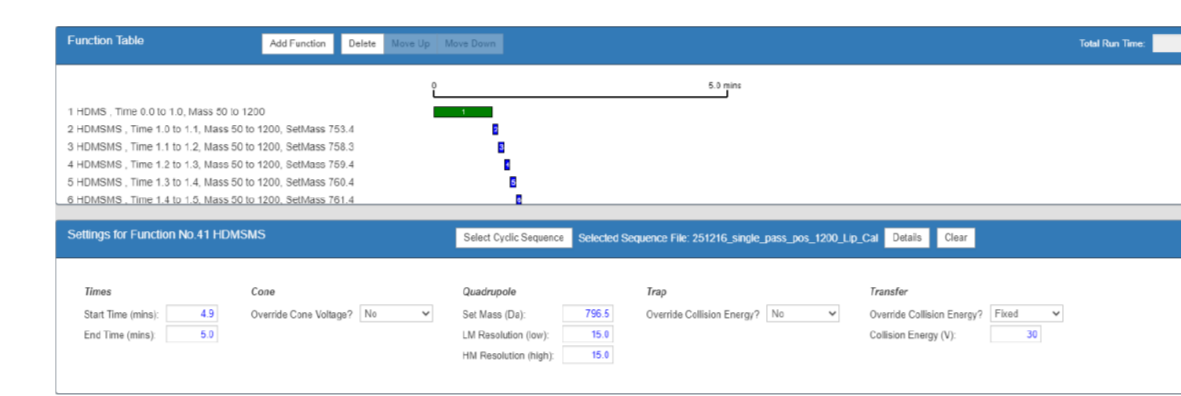


Figure 4. Proof of principle method design for collecting cIM-MS and Q-cIM-MS/MS

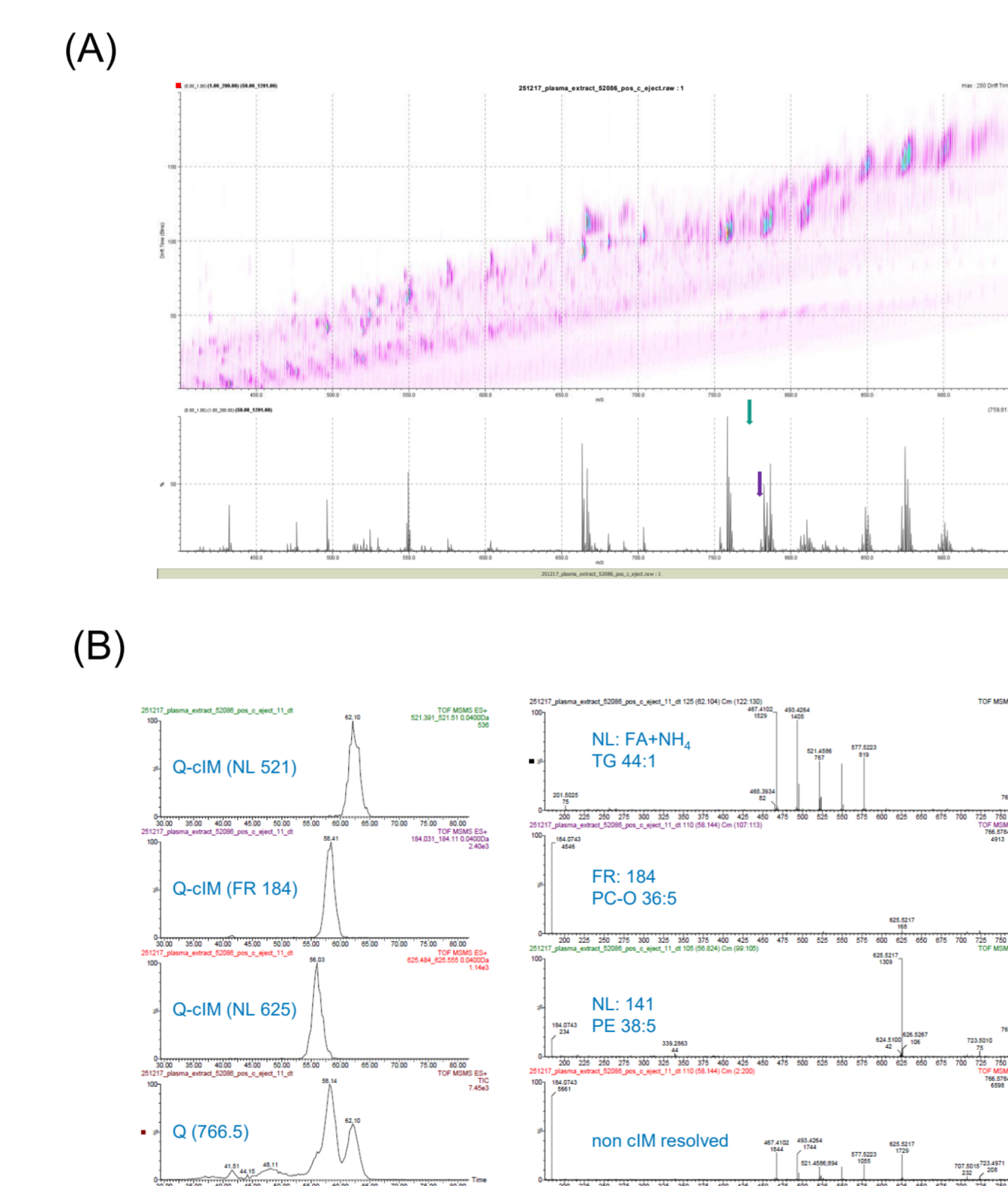


Figure 5. (A) Lipid plasma extract m/z vs. drift time cIM-MS profile, and (B) fragment ATDs of m/z 766.5 (left) and non- and cIM-resolved MS/MS spectra (right), including lipid annotation.

Ion mobility–resolved association of Q-cIM-MS/MS and low-energy cIM-MS enables minimization of signal overlap from isobaric and isomeric lipid species in a shotgun lipidomics approach. The data presented here are based on single-pass experiments.

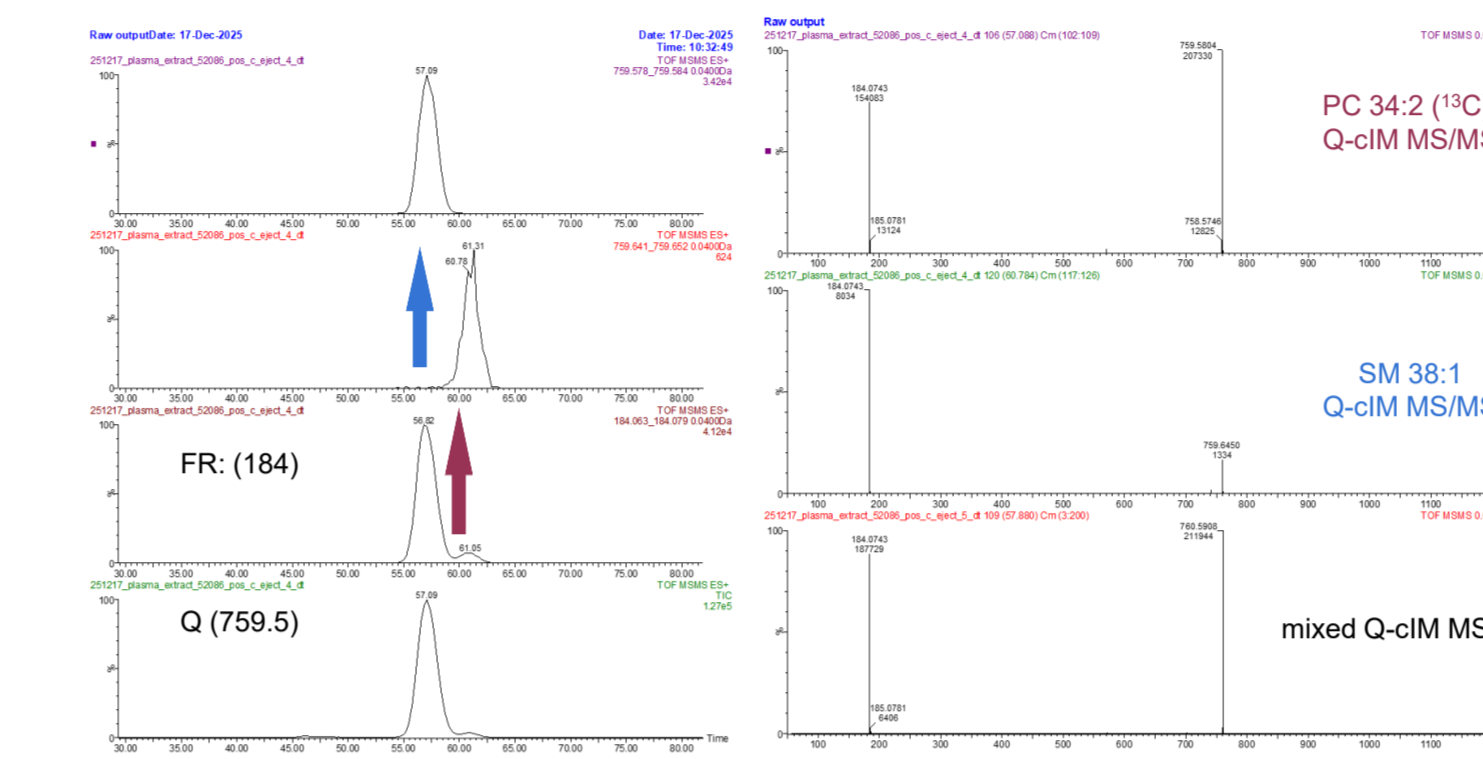


Figure 6. Non-resolved and cIM resolved ATDs of m/z 759.5 (left) and the corresponding MS/MS spectra (right), showing that isomeric overlap of FR: 184.07 can be IM resolved.

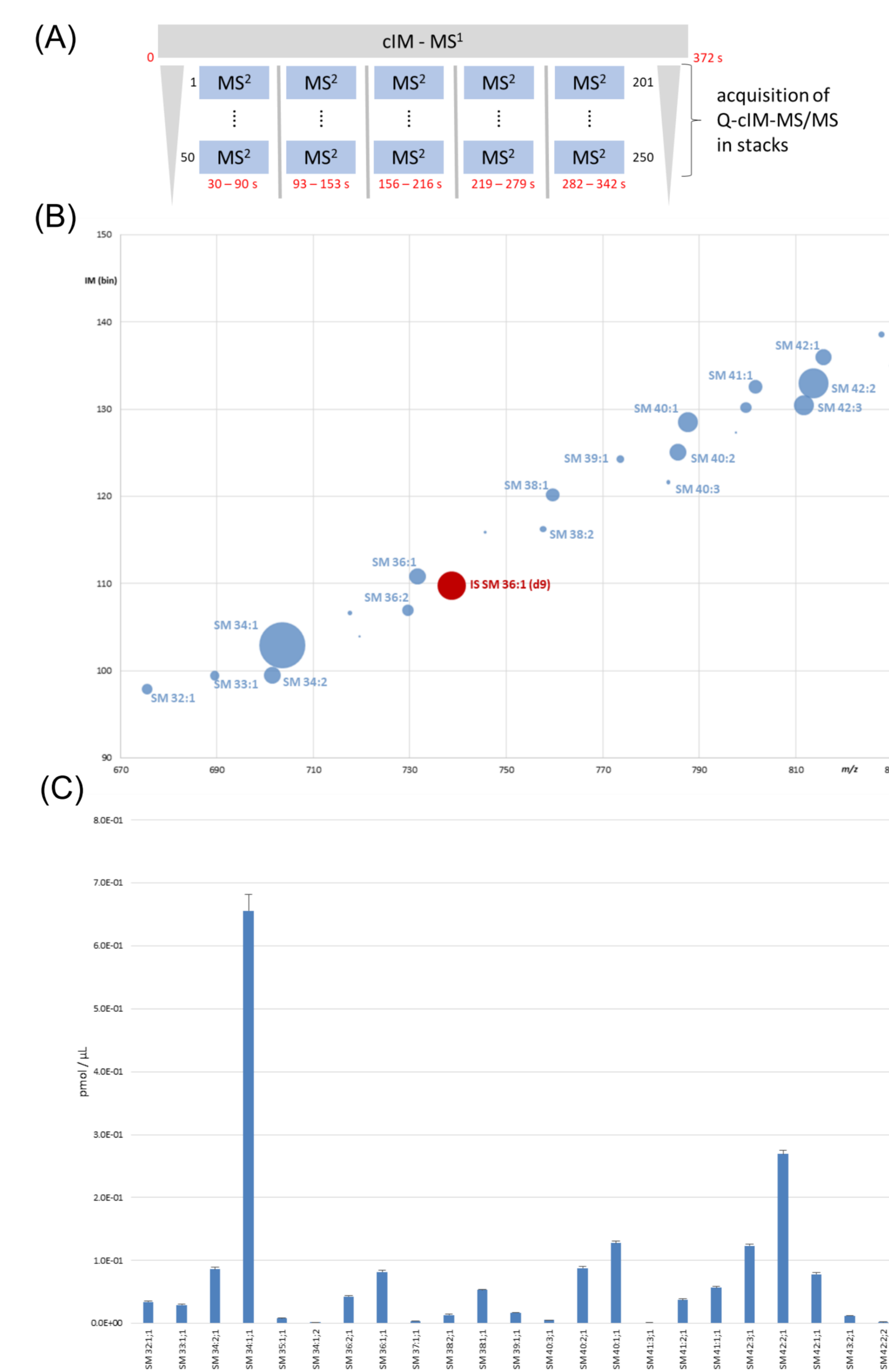


Figure 7. (A) Design of a Q-cIM-MS/MS direct infusion experiment, in which 250 Q-cIM-MS/MS were acquired within 6 min. (B) and (C) IM confirmed SM species.

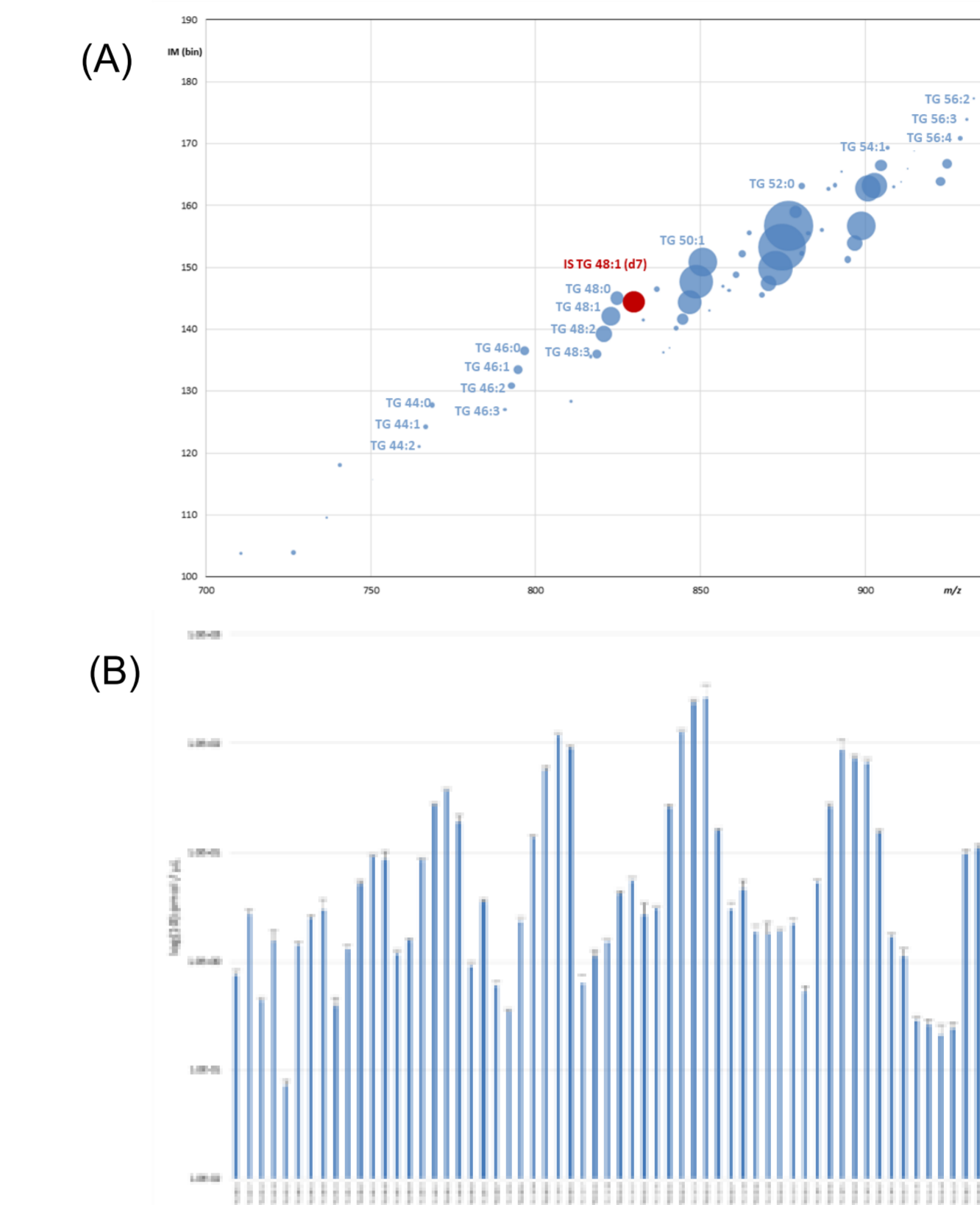


Figure 8. (A) TG species in 2D representation (m/z , IM). (B) Quantitation results.


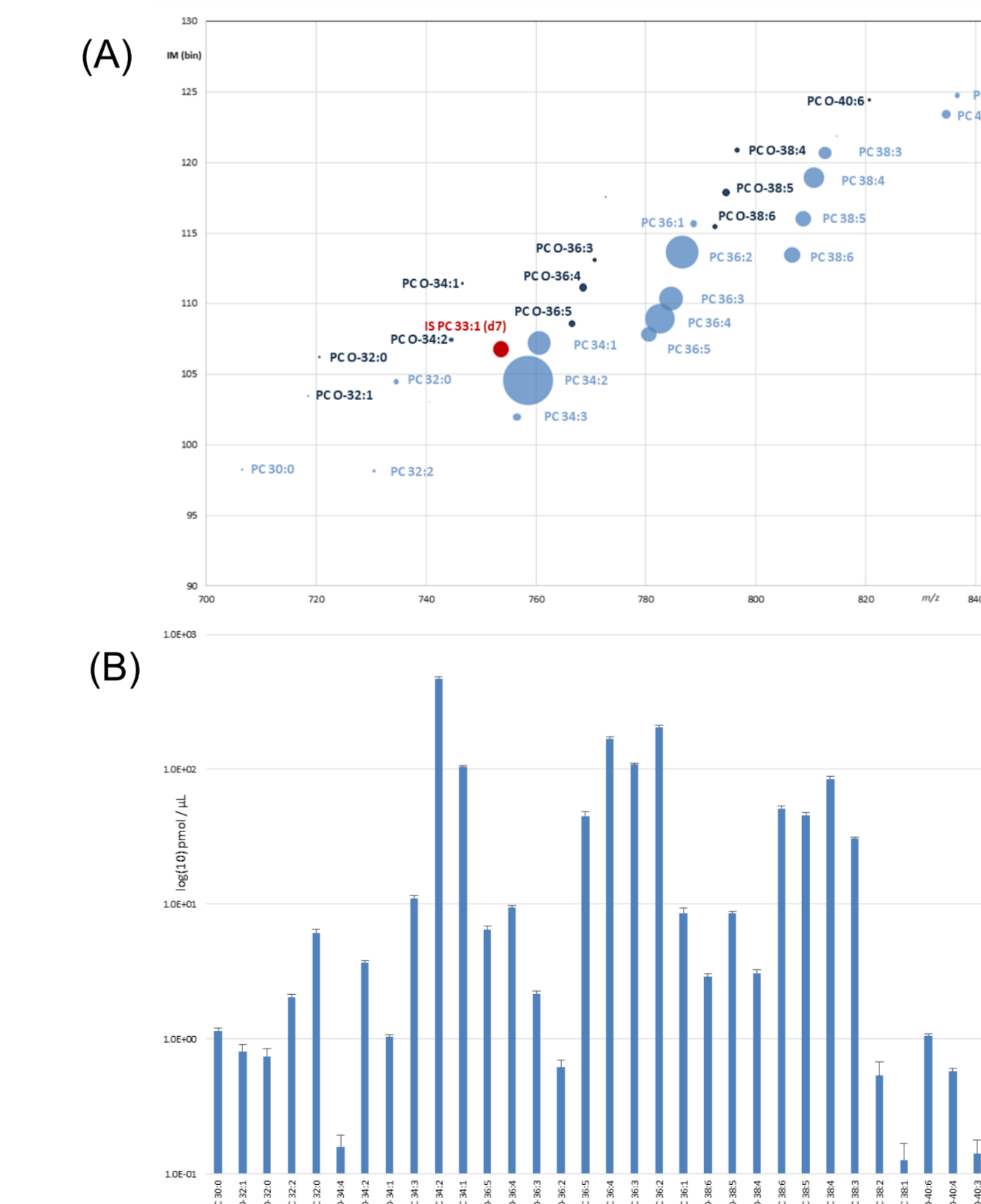


Figure 9. (A) Phosphatidylcholine (PC) and ether-linked phosphatidylcholine (PC-O) are separated in 2D. (B) Overall PC lipid species profile.

Conclusions

- Q-cIM-MS/MS acquisitions produces MS/MS spectra that resolve complex isomeric and isobaric overlaps
- Multidimensional peak detection was performed as part of the data processing pipeline used for both method creation and lipid annotation.
- The presented workflow establishes a framework that can be extended to larger sample cohorts
- A total of 166 lipids from eight classes were quantified using positive-ion shotgun lipidomics in a proof-of-principle experiment

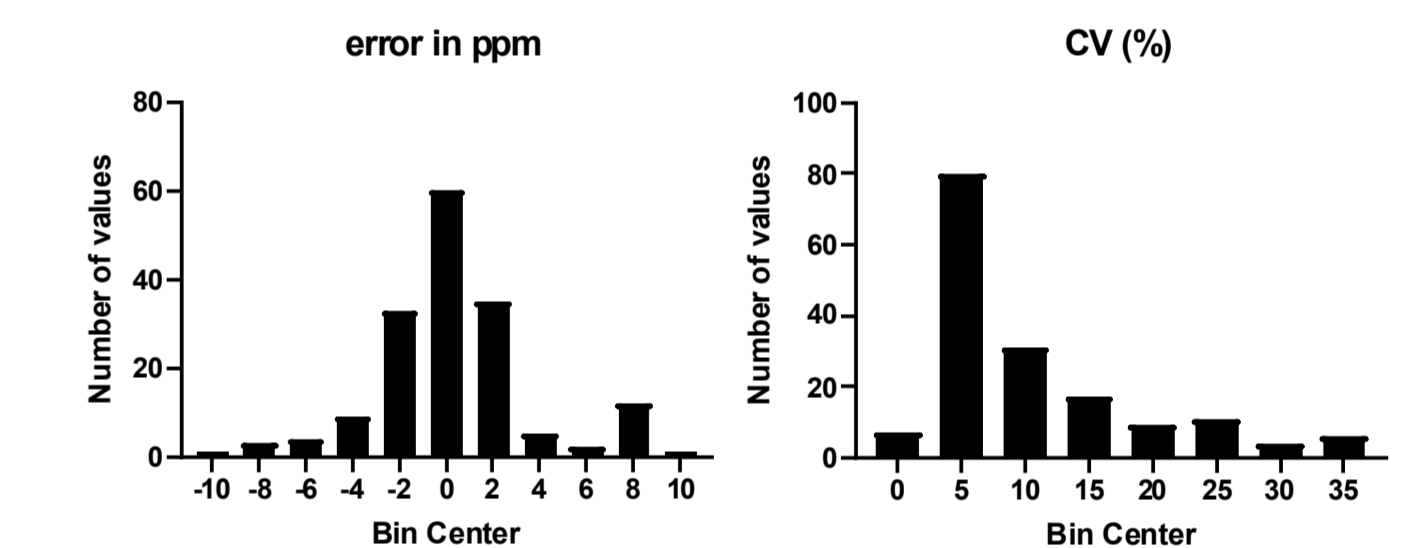

Outlook

- Integration of Q-cIM-MS/MS in positive and negative ion modes
- Increased Q-cIM-MS/MS coverage via inclusion lists
- Automated generation of 2D feature lists
- LipidXplorer updates for multidimensional MasterScan generation and lipid identification

Acknowledgement

Funding for this research in the DS Lab was supported by the German Network for Bioinformatics Infrastructure (de.NBI)

References

1. Giles et al. A Cyclic Ion Mobility–Mass Spectrometry System. *Anal Chem.* 2019
2. Matyash et al. Lipid extraction by methyl-tert-butyl ether for high-throughput lipidomics. *JLR* 2008
3. Herzog et al. A novel informatics concept for high-throughput shotgun lipidomics based on the molecular fragmentation query language. *Genome Biol.* 2011
4. <https://lifs-tools.org/tools/ixpostman.html>

Python is a registered trademark of the Python Software Foundation. SELECT SERIES, Cyclic, MassLynx, DriftScope and Waters are trademarks of Waters Corporation or its affiliates. Triversa and NanoMate are trademarks of Advion Interchim Scientific