



## INTRODUCTION

- Native top-down protein analysis is advancing the characterisation of proteoforms in a biological context
- Spectral resolution is often limited by incomplete desolvation of ions from endogenous matrices, salts, and solubilizing detergents<sup>1,2</sup>
- High quality tandem mass spectra remain challenging to obtain for low abundance proteoforms and large oligomeric assemblies<sup>3</sup>
- Here, a declustering ion guide improves protein resolution, and Wideband Enhancement (WBE) boosts tandem MS sensitivity
- NISTmAb was used to develop methods and optimize declustering, and the declustering guide is shown to enable pseudo-MS<sup>3</sup> experiments for top-down sequencing of streptavidin.
- Furthermore, we show the guide to facilitate the release of bacteriorhodopsin from a detergent micelle to enable characterization of bound lipids

## METHODOLOGY

### Samples

- NISTmAb standard (Pierce)
- Streptavidin (Pierce, 21125)
- Bacteriorhodopsin (bR) from *halobacterium salinarum* (Sigma-Aldrich)

Samples were electrosprayed from 200 mM ammonium acetate using 2 μm internal diameter borosilicate glass nanocapillaries (WPI LLC, FL, USA). To prepare native bacteriorhodopsin, lyophilized samples were resuspended, vortexed and buffer exchanged in 40 mM β-D-glucopyranoside (OG) detergent (Sigma-Aldrich) and 200 mM ammonium acetate.

### Mass Spectrometry

All experiments were performed with a Waters™ Cyclic™ IMS P20 Mass Spectrometer operating in positive nESI mode, equipped with a Dynamic Field Declustering (DFD) device (Figure 1). Data were acquired in MassLynx™ v4.2 Software, processed using the top-down workflow in waters\_connect™ UNIFI™ software.

## REFERENCES

- Cheng D, Guo Y, Lyu J, et al. Advances and challenges in preparing membrane proteins for native mass spectrometry. *Biotechnology Advances*. 2025;78:108483.
- Tamara S, den Boer MA, Heck AJR. High-Resolution Native Mass Spectrometry. *Chem Rev*. 2022;122(8):7269-7326.
- Hoang C, Uritboonthai W, Hoang L, et al. Tandem Mass Spectrometry across Platforms. *Anal Chem*. 2024;96(14):5478-5488.
- Inada M, Kinoshita M, Matsumori N. Archaeal Glycolipid S-TGA-1 Is Crucial for Trimer Formation and Photocycle Activity of Bacteriorhodopsin. *ACS Chem Biol*. 2020;15(1):197-204.

## ACKNOWLEDGEMENTS & STATEMENTS


**Conflict of Interest statement:** All authors are employees of Waters Corporation who manufacture and sell the products highlighted in this work.

Waters, Cyclic, MassLynx, waters\_connect and UNIFI are trademarks of Waters Corporation or its affiliates.

The authors are grateful to the British Mass Spectrometry Society, American Society of Mass Spectrometry and Waters Corporation for providing financial support and grants to present this work.

## TECHNOLOGY


### Dynamic Field Declustering

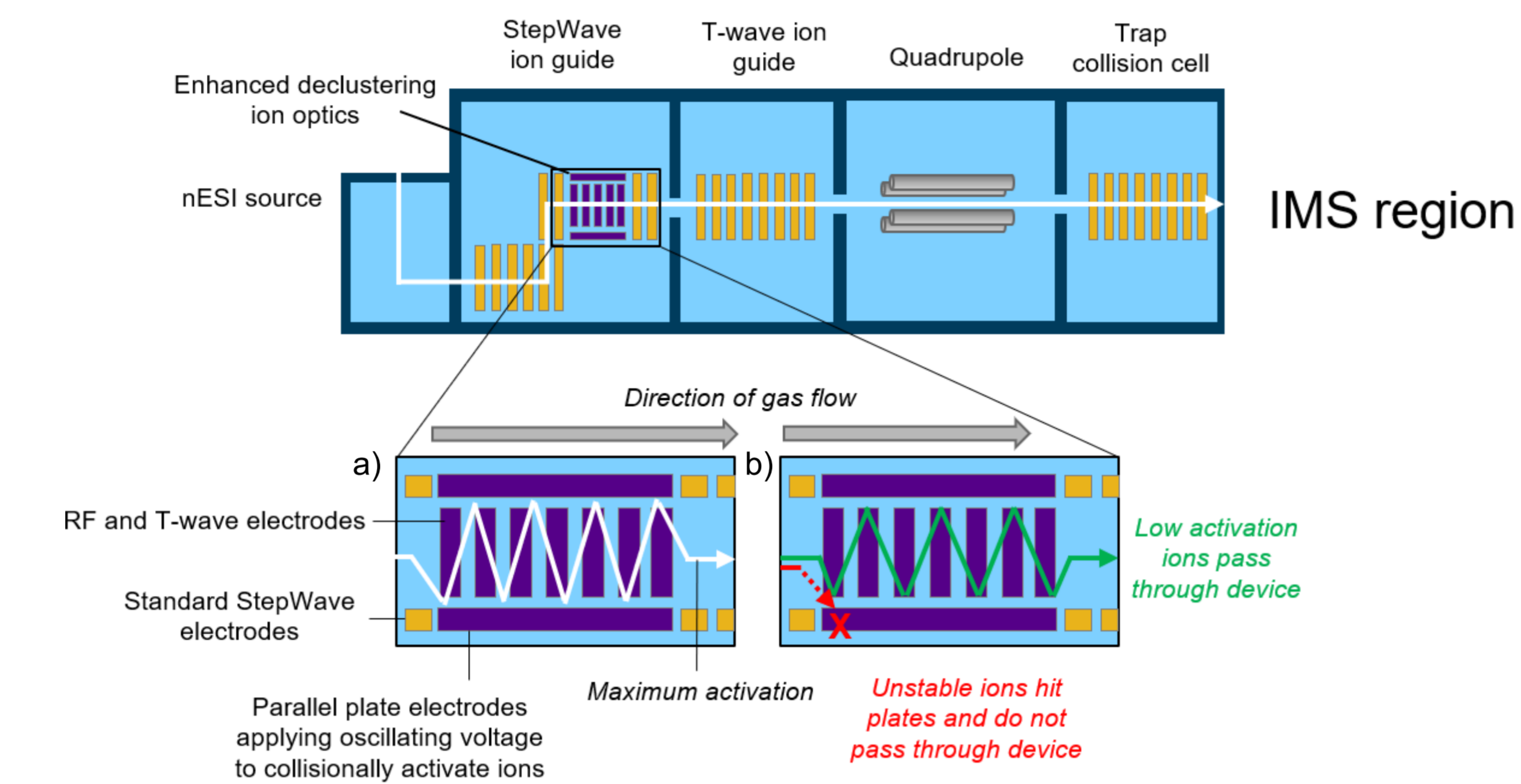

**Figure 1.** Schematic of the DFD device. Pre-IMS region is shown (TOP), highlighting the location of the declustering ion guide. **a)** Demonstrating ion movement with maximum activation, whereby the longest distance is travelled. **b)** Highlighting how ions are separated by different paths: unstable pathway, and low activation pathway, which continue into the quadrupole and beyond for detection.

### Wideband Enhancement


In Time of Flight (TOF) operation, incomplete transmission from a continuous beam limits sensitivity; WBE increases duty cycle by synchronising ion release with mobility-separated ion packets, improving sampling efficiency across a broad mass range.

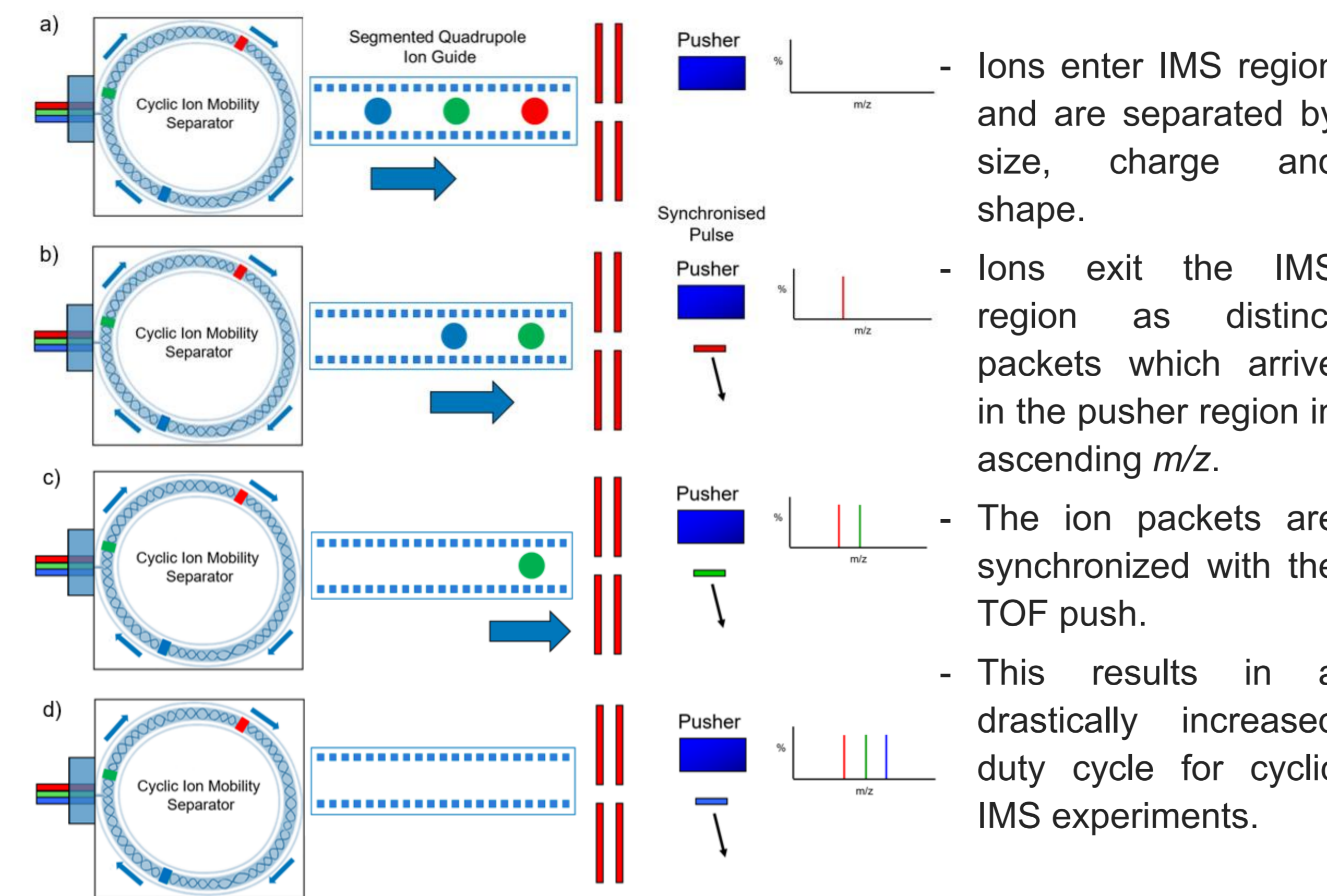

**Figure 2.** Schematic of WBE mode of operation and the corresponding spectra produced


## RESULTS

### Optimizing declustering with NISTmAb

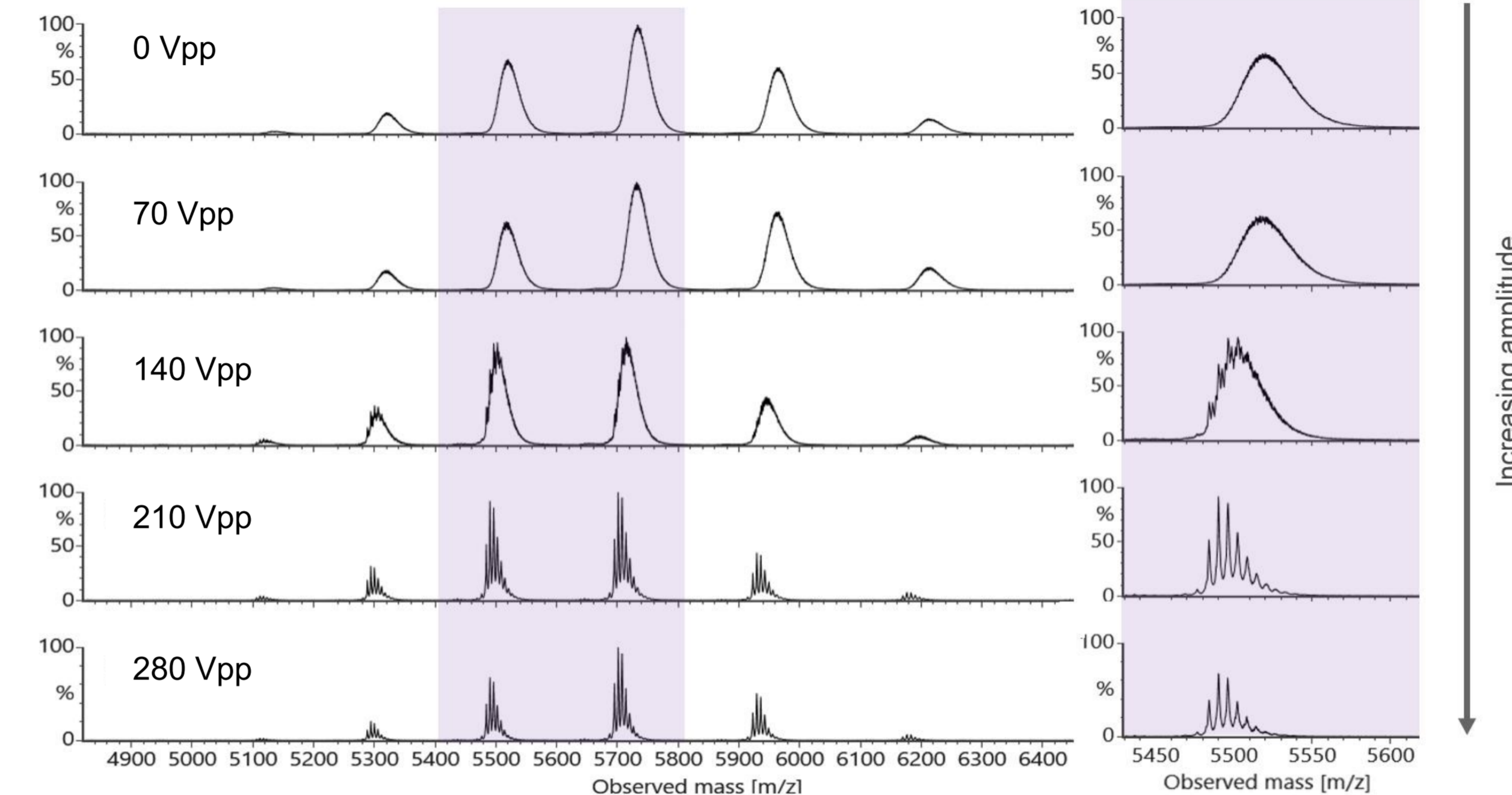


**Figure 3.** Nano-electrospray ionization spectra for NISTmAb antibody at increasing DFD amplitudes and fixed frequency (60 kHz). Increasing amplitudes from 0-280 Vpp improved declustering and apparent peak resolution of NISTmAb glycoforms, demonstrating that DFD can simply and efficiently desolvate and decluster native samples with minimal tuning required.

### Top-Down MS of Streptavidin

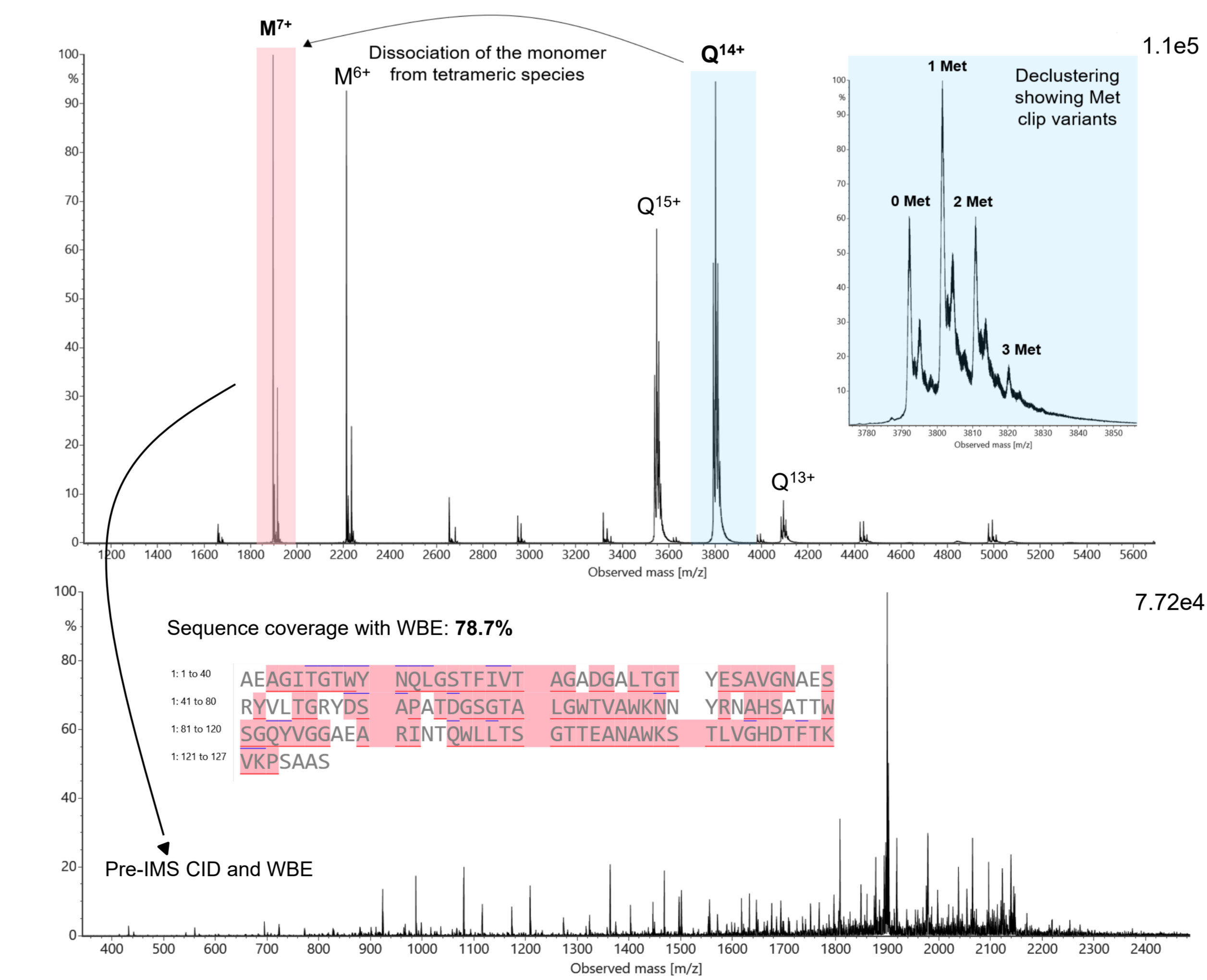


**Figure 4.** Top-down MS of native streptavidin. (TOP) DFD set at an amplitude of 175 Vpp and a 60 kHz frequency declustered the tetramer (Q), improving apparent peak resolution to show the Met clip variants (blue), in addition to partially dissociating the 7+ monomer (M) (red). (BOTTOM) The monomer was quadrupole selected and fragmented by CID in a pseudo-MS<sup>3</sup> experiment for top-down analysis, using WBE to produce a high-quality spectrum, achieving 78.7% coverage.

### Native bacteriorhodopsin

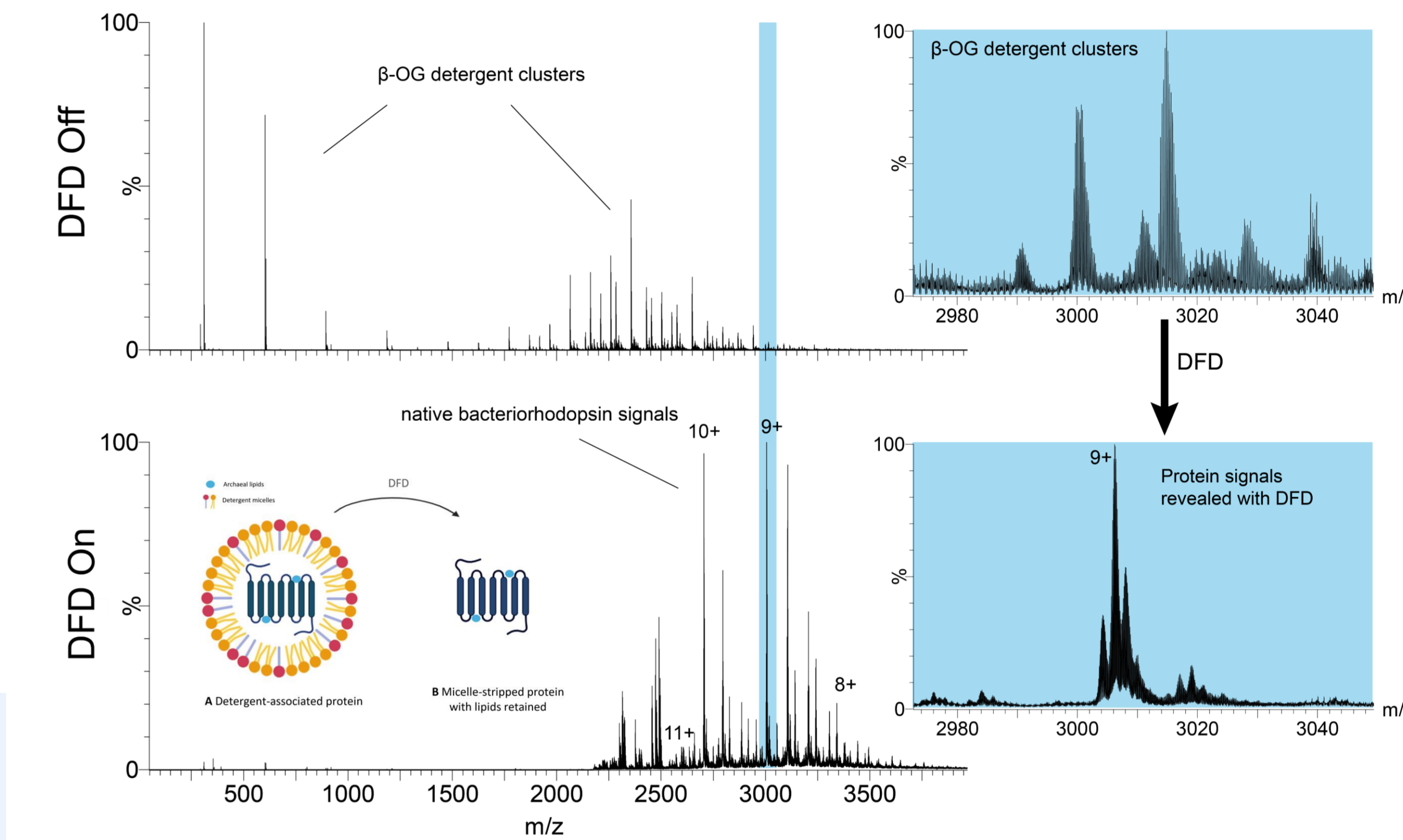

**Figure 5.** Utilizing DFD to study native bR releases the protein from the micelle to reveal protein signals on the native mass spectra. The blue inset highlights the 9+ charge state of monomeric native bR, demonstrating how DFD effectively declusters the protein signals from detergent to reveal the isotopically resolved monomer.

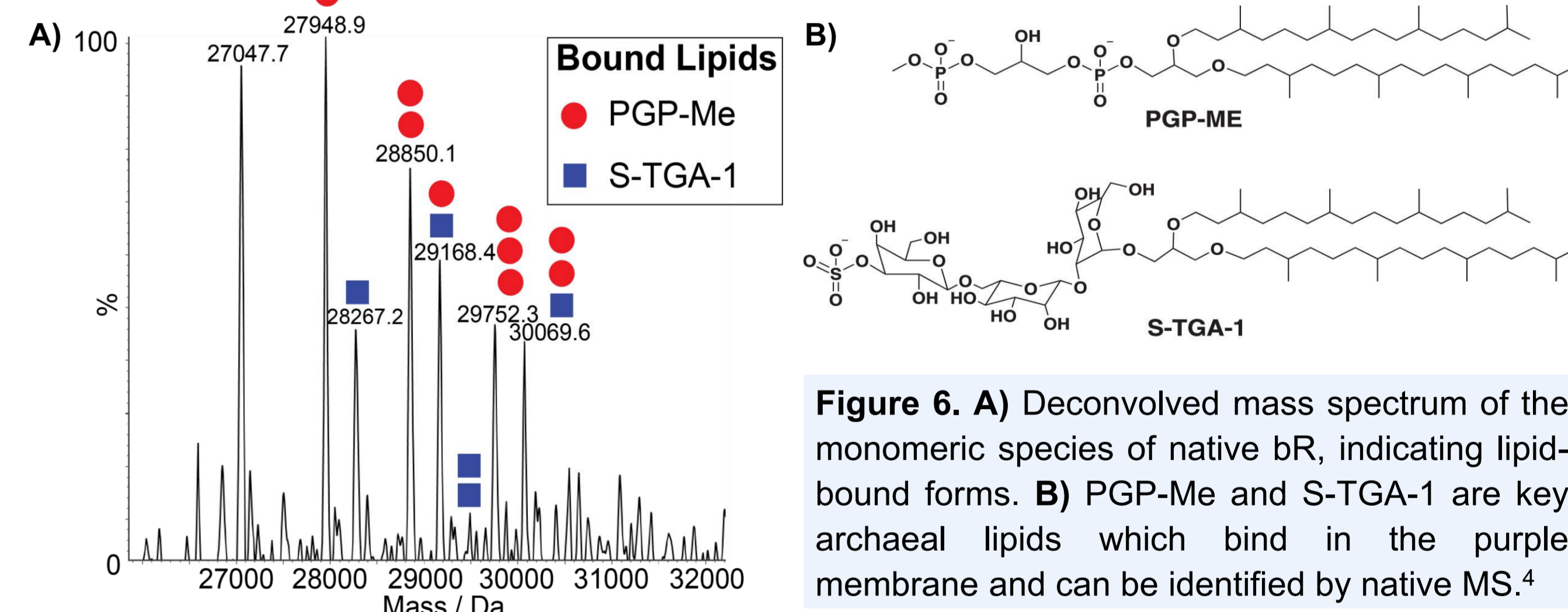

**Figure 6.** **A)** Deconvolved mass spectrum of the monomeric species of native bR, indicating lipid-bound forms. **B)** PGP-Me and S-TGA-1 are key archaeal lipids which bind in the purple membrane and can be identified by native MS.<sup>4</sup>

## CONCLUSIONS & FUTURE WORK

### Conclusions:

- DFD streamlines native MS by efficiently removing salt/solvent adducts and by releasing membrane proteins from detergent micelles.
- WBE increases sensitivity to deliver high quality top-down data in less time.
- In combination with DFD, WBE enables pseudo-MS<sup>3</sup> workflows, yielding sequence level information directly from native complexes.

### Future work:

- This work establishes an enhanced analytical framework for native proteins, providing a robust platform for continued investigation.
- Ongoing studies focus on bacteriorhodopsin, with development of sample preparation strategies to support trimer detection and enable top-down analysis of the native complex