

# Analysis of Aliphatic Alcohols by Ligand-Exchange Chromatography

## Application Note

Chemical

### Author

Stephen Ball  
Agilent Technologies, Inc.

### Introduction

This application note demonstrates how an Agilent Hi-Plex H column can be used to separate aliphatic alcohols.



**Agilent Technologies**

## Materials and Reagents

Column	Agilent Hi-Plex H (8% crosslinked), 7.7 × 300 mm, 8 μm (p/n PL1170-6830)
Mobile phase	100% DI H <sub>2</sub> O
Flow rate	0.6 mL/min
Temperature	40 °C
Detector	RI

## Conclusion

Using only pure HPLC-grade water as eluent, the Agilent Hi-Plex H column is capable of separating a range of aliphatic alcohols. In addition to those shown in Figure 1, it may also be possible to separate a much wider range of this type of compound. Molecular weight and degree of branching are critical factors in determining the amount of retention on a Hi-Plex H column.

## For More Information

These data represent typical results. For more information on our products and services, visit our Web site at [www.agilent.com/chem](http://www.agilent.com/chem).

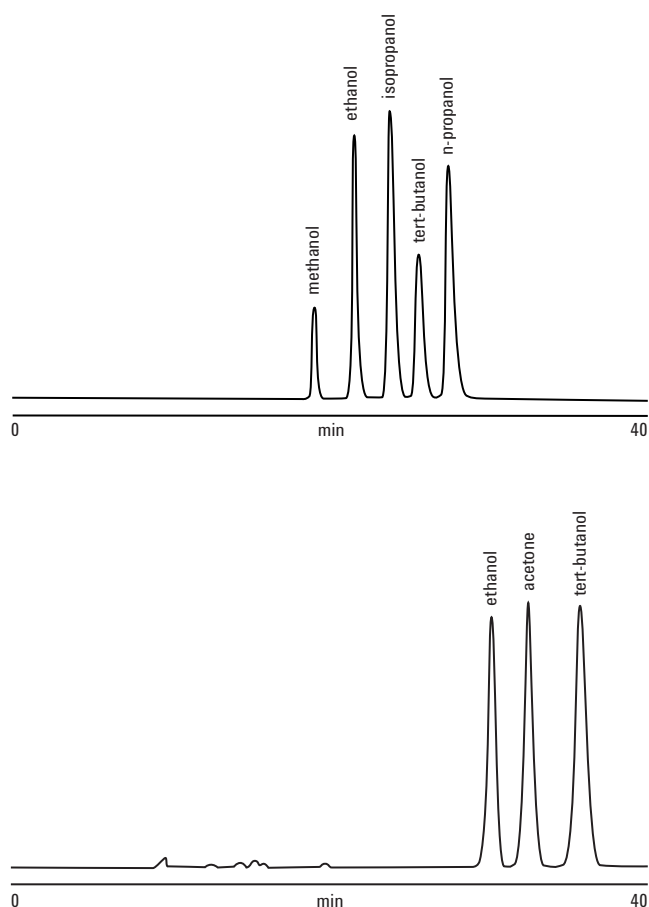


Figure 1. Separation of different aliphatic compounds on an Agilent Hi-Plex H column.

[www.agilent.com/chem](http://www.agilent.com/chem)

Agilent shall not be liable for errors contained herein or for incidental or consequential damages in connection with the furnishing, performance, or use of this material.

Information, descriptions, and specifications in this publication are subject to change without notice. This publication was originally published in 2009.

© Agilent Technologies, Inc., 2011  
Published in USA, June 30, 2011  
SI-02030



**Agilent Technologies**