

## Application News

Ultra High Performance Liquid Chromatograph Nexera™ XS inert

# Improvement of Oligonucleotide Peak Shape Using Automatic Pretreatment Function (Co-injection)

Shinichi Fujisaki, Momoka Hayashida

### User Benefits

- ◆ Sample dilution can be automated using the automatic pretreatment function “co-injection” for purified fractions of oligonucleotides containing high concentrations of salts.
- ◆ The automatic pretreatment function “co-injection” enables the simultaneous injection of a sample and desired solvents.

### ■ Introduction

In HPLC analysis, the composition of the sample solvent is critical for achieving proper peak shape. If the sample solvent is a stronger eluting solvent than the mobile phase, sample band condensation at the column inlet may be insufficient, leading to sample band broadening. For instance, increasing the organic solvent ratio in the sample solvent to dissolve low-polar compounds may deteriorate the peak shapes of early-eluting compounds in reversed-phase chromatography. Oligonucleotides are synthesized using the phosphoramidite method and subsequently purified by anion exchange chromatography or reversed-phase ion-pair chromatography (RP-IP). Purified fractions from anion exchange chromatography often contain high concentrations of salts, such as sodium chloride (NaCl) and sodium bromide (NaBr). When these purified fractions are analyzed using RP-IP, the high salt concentration in the sample solvent may interfere with the ion-pair formation capacity of the oligonucleotides. This interference can result in improper peak retention and adversely affect the chromatogram. This article introduces a case study demonstrating an improvement in oligonucleotide peak shape by suppressing the sample solvent effect through simultaneous injection with a desired solvent. This automatic pretreatment function “co-injection” is a standard feature included in the autosamplers of Nexera series.

### ■ “Co-injection” Function

The co-injection function, equipped in autosamplers, provides automatic pretreatment. A specified amount of reagent or solvent from any vial can be mixed with the sample solution in the needle and injected (Fig. 1).

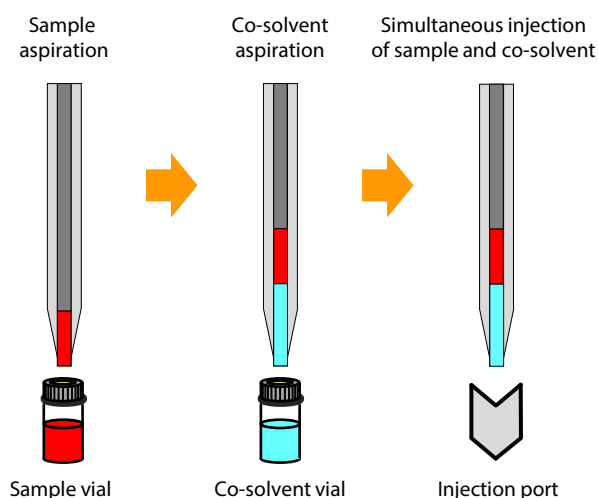


Fig. 1 Flow Diagram of Co-injection Sequence

### ■ Analytical Conditions

The analytical conditions are shown in Table 1. Two types of salts, 3 mol/L NaCl and 3 mol/L NaBr, were added to the sample solvent to simulate purified fractions containing high concentrations of salts. To assess the effects of co-solvents and co-injection volumes on peak shape, four different co-solvents were used: water and three solutions containing HA (hexylamine) at concentrations of 10, 30, and 50 mmol/L, each prepared in 100 mmol/L HFIP (1,1,1,3,3,3-hexafluoro-2-propanol). Additionally, three injection volumes (1, 3, and 5  $\mu$ L) were evaluated. The target sample was one of the models of oligonucleotide therapeutics synthesized in the AMED project<sup>\*1</sup>.

\*1 Grant Number: JP21ae0121022, JP21ae0121023, JP21ae0121024

Table 1 Analytical Conditions

System	: Nexera XS inert (Method Scouting System)
Column	: Accura Triart Bio C18 <sup>*2</sup> (100 mm $\times$ 2.1 mm I.D., 1.9 $\mu$ m)
Sample solvents	: 20 mM Tris-HCl (pH 8.5) containing 3 mol/L NaCl : 20 mM Tris-HCl (pH 8.5) containing 3 mol/L NaBr
Co-solvents	: Water : 100 mmol/L HFIP <sup>*3</sup> and 10 mmol/L HA <sup>*4</sup> in water : 100 mmol/L HFIP and 30 mmol/L HA in water : 100 mmol/L HFIP and 50 mmol/L HA in water
Co-injection volumes	: 1, 3, 5 $\mu$ L
Mobile phases	
Pump A	: 100 mmol/L HFIP and 10 mmol/L HA in water
Pump B	: 100 mmol/L HFIP and 10 mmol/L HA in Methanol
Temperature	: 60 °C
Injection volume	: 1 $\mu$ L
Flow rate	: 0.2 mL/min
Time program (%B)	: 25 % (0-2 min) $\rightarrow$ 67 % (26 min) $\rightarrow$ 100 % (26-27 min) $\rightarrow$ 25 % (27-35 min)
Detection	: 260 nm (SPD-M40, UHPLC inert cell)

\*2 YMC.CO., LTD.

\*3 1,1,1,3,3,3-hexafluoro-2-propanol

\*4 Hexylamine

### ■ Improved Peak Shape Using Co-injection Function

Chromatograms obtained without the co-injection function are shown in Fig. 2(2) (sample solvent: NaCl) and Fig. 4(2) (sample solvent: NaBr), respectively. The ion-pair formation ability was suppressed due to the high concentration of salts in the sample solvents. Consequently, the oligonucleotide was not properly retained, and part of the peak was eluted as a split peak. On the other hand, simultaneous co-injection of pure water resulted in increased peak areas for the full-length product (FLP) (Fig. 2 (3)–(5) and Fig. 4 (3)–(5)), confirming that it effectively provided appropriate peak retention. When the co-injection volume was 3  $\mu$ L or more (Fig. 2(4), (5) and Fig. 4(4), (5)), split peaks were no longer observed. The relationship between the co-solvents, co-injection volumes, and the normalized peak areas (%) of FLP is shown in Fig. 3 (sample solvent: NaCl) and Fig. 5 (sample solvent: NaBr), respectively. Regardless of the co-injected solvent type, the normalized peak areas of FLP remained nearly constant when the co-injection volume was 3  $\mu$ L or more. This suggests that co-injection is effective in suppressing the peak splitting phenomenon caused by the sample solvent effect.

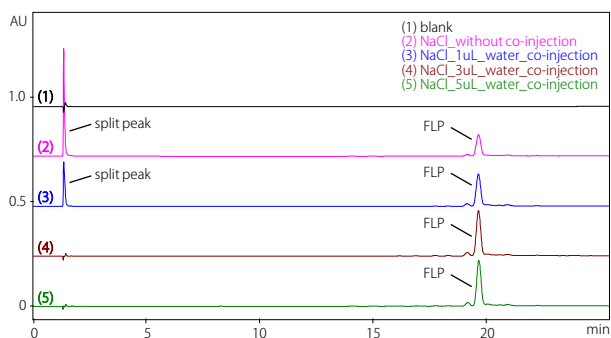


Fig. 2 Comparison of Chromatograms with/without Co-injection (sample solvent : NaCl)

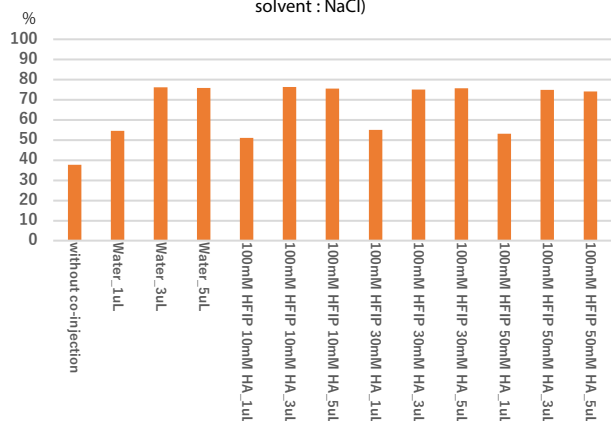


Fig. 3 Normalized FLP Peak Areas (sample solvent : NaCl)

\* normalization based on sum of all peaks including FLP related impurities

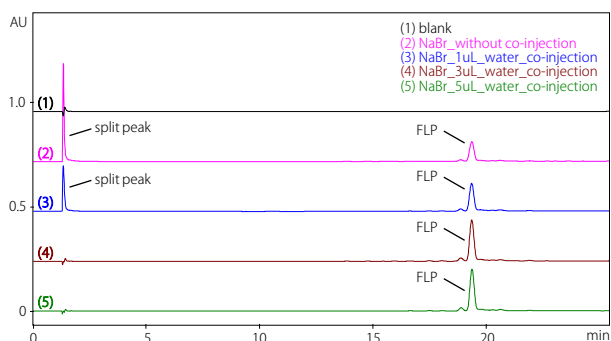


Fig. 4 Comparison of Chromatograms with/without Co-injection (sample solvent : NaBr)

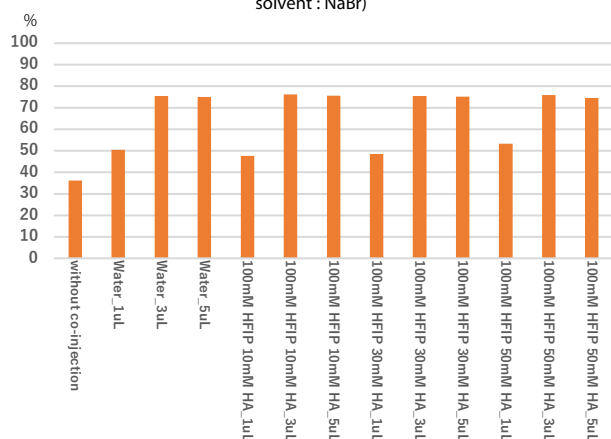


Fig. 5 Normalized FLP Peak Areas (sample solvent : NaBr)

\* normalization based on sum of all peaks including FLP related impurities

LabSolutions, Nexera, DUIS, and Shim-pack Scepter are trademarks of Shimadzu Corporation or its affiliated companies in Japan and/or other countries.

## Conclusion

When purified fractions of oligonucleotides containing high concentrations of salt are analyzed by RP-IP, the ion-pair formation capacity is suppressed, and peaks may not be properly retained, resulting in partial early elution as a split peak. Pretreatment, such as desalination or dilution of samples, is usually required before analysis, but it is time-consuming. The co-injection function, equipped in Nexera series autosamplers, allows for the simultaneous injection of a desired solvent with purified fractions, providing automatic pretreatment that enables the direct analysis of oligonucleotides in purified fractions with high salt concentrations. In particular, when many purified fractions need to be analyzed, significant labor savings can be achieved because the necessary pretreatment procedures are automated for all fractions. The optimal co-injection solvent and volume depend on the analytical conditions, such as oligonucleotide sequence and mobile phase composition. Therefore, optimization of injection conditions is necessary on each case, and the co-injection function enables more efficient comparative investigations under multiple different injection conditions.

## Acknowledgments

A part of this work was supported by AMED under Grant Number JP21ae0121022, JP21ae0121023, JP21ae0121024 (Project leader: Satoshi Obika).



# SHIMADZU

Shimadzu Corporation

www.shimadzu.com/an/

For Research Use Only. Not for use in diagnostic procedures.

This publication may contain references to products that are not available in your country. Please contact us to check the availability of these products in your country.

The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. See <http://www.shimadzu.com/about/trademarks/index.html> for details.

Third party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

Shimadzu disclaims any proprietary interest in trademarks and trade names other than its own.

The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.

01-00848-EN

First Edition: May, 2025

› Please fill out the survey

## Related Products

Some products may be updated to newer models.



› Nexera XS inert  
Ultra High Performance Liquid  
Chromatograph

## Related Solutions

› Pharmaceutical and  
Biopharmaceutical

› Oligonucleotide and  
mRNA Therapeutics

› Price Inquiry

› Product Inquiry

› Technical Service /  
Support Inquiry

› Other Inquiry