



Batch copy or export of
multiple images from MS
Image List

Export all displayed images → Right-click on "Outside images".

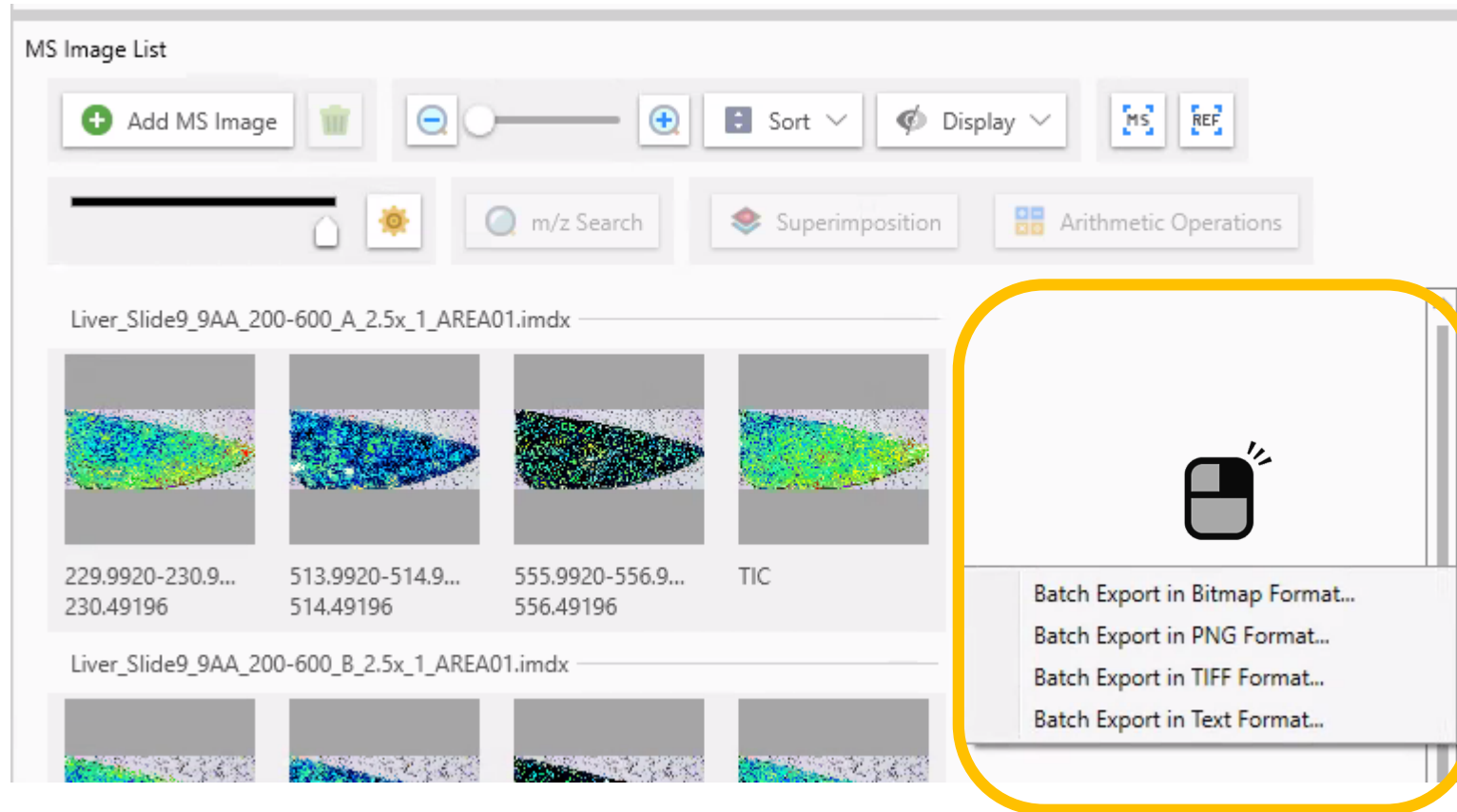

※In this case, the "MS Image Setting" will not be applied.


Copy or export the selected images →
Right-click “on the image”.

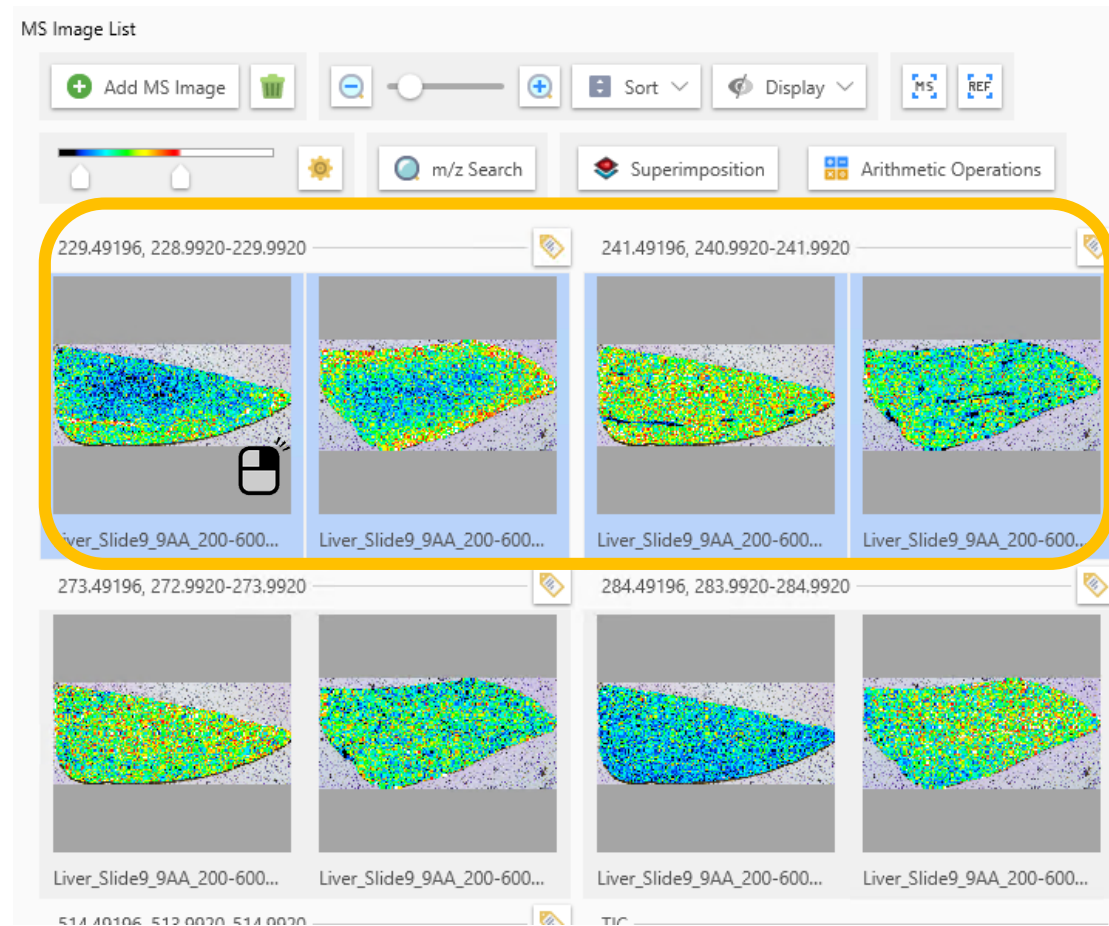

Select multiple images


Select from the context menu

The screenshot displays the 'MS Image List' software interface. At the top, there is a toolbar with various icons for adding, deleting, zooming, sorting, and displaying images. Below the toolbar, there are several control panels including a color scale, a search bar for 'm/z', and buttons for 'Superimposition' and 'Arithmetic Operations'. The main area shows a grid of mass spectrometry images. A context menu is open over one of the images, listing the following options: 'Delete', 'Batch Copy Sort Group', 'Batch Export Sort Group...', 'Copy Selected MS images', and 'Export Selected MS images...'. The images in the grid are labeled with m/z ranges and file names, such as 'Liver_Slide9_9AA_200-600...'.

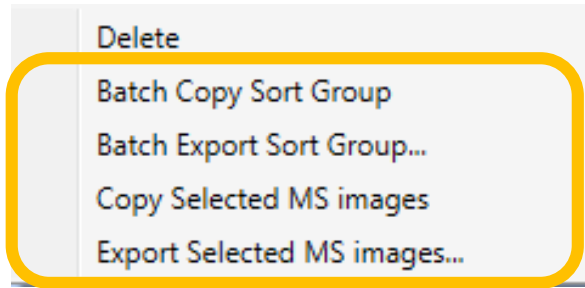

- “Batch Copy Sort Groups”
 - Copies all images in the group that contains the right-clicked image.
 - *The group method can be changed in "Sort".
- “Batch Export Sorted Group”
 - Batch export the images of the group that contains the right-clicked image.
- “Copy selected MS images”
- “Export selected MS images”

Change output format

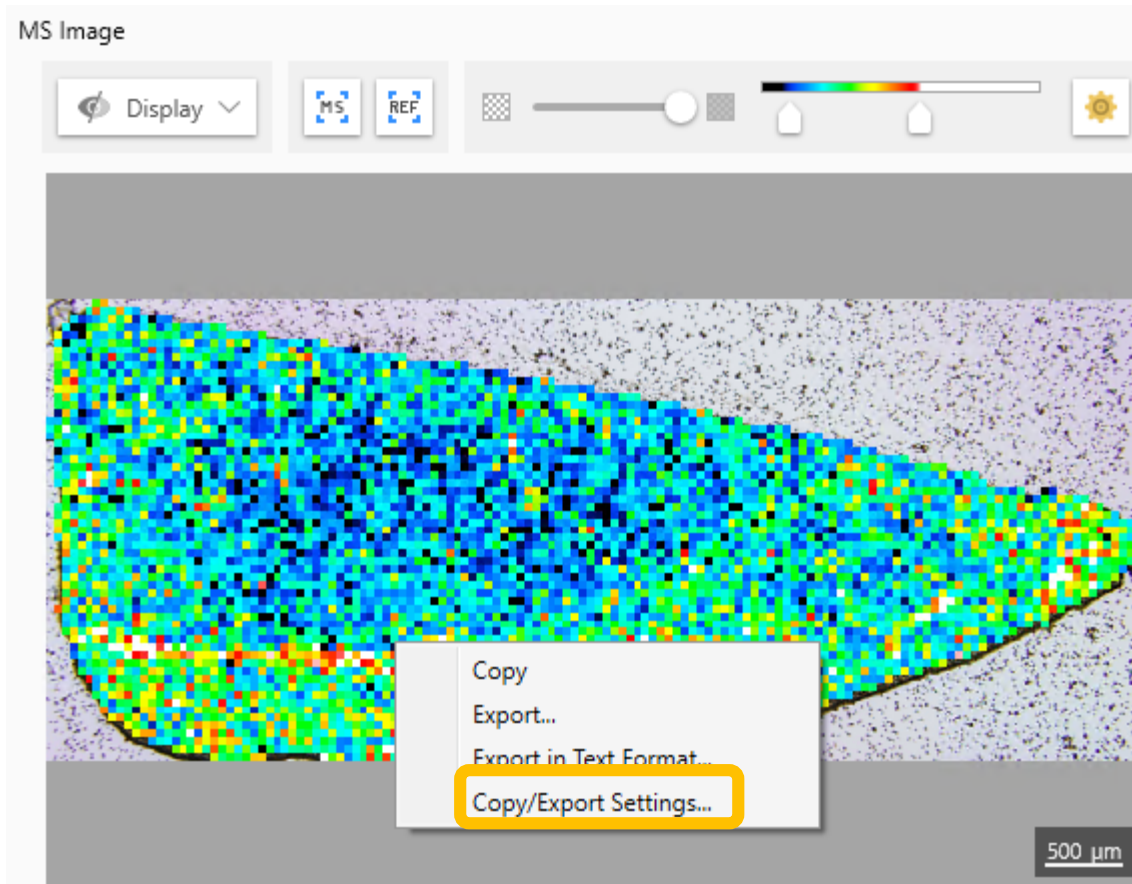

The output format can be changed from "Copy/Export Settings" under "MS Image".

