CoreFocus

Syringe Filter

OPERATION GUIDE

Introduction

Thank you very much for purchasing this product. This disposable, non-sterile syringe filter is designed and manufactured to efficiently remove particulate impurities from sample solutions. It is suitable for pretreatment of samples for analytical instruments such as Ion Chromatography and HPLC, as well as for clarification or particle removal of various other liquid samples. A wide range of types is available to meet the volume and properties of your samples. This guide contains important information required for proper use of this product. Please read it carefully before use.

Handling Precautions

This product is for research use only and is not a medical device.

- The product is non-sterile. Do not use it for sterile filtration.
- To prevent contamination, do not filter more than one type of sample with a single filter unit.

Specifications

Filter Size	4 mm	13 mm	25 mm
Pore size (μm)	0.2 / 0.45		
Material	Olefin-based polymer		
Effective filtration area (cm ²)	0.07	0.8	4
Max. operating pressure (MPa)	0.49		
Max. operating temperature (°C)	100		
Hold-up volume	≤10 μL	≤30 μL	≤100 μL
Inlet connector	Luer lock		
Outlet connector	Luer slip		
Quantity	100 pieces		

Sample volume guideline is as follows.

Filter Size	Sample Volume	Recommended Syringe Volume
4 mm	≤1 mL	1 mL, 2 mL
13 mm	1 - 10 mL	2 mL, 5 mL, 10 mL
25 mm	10 - 50 mL	10 mL, 20 mL, 50 mL

Operating Precautions

- Use a syringe with capacity sufficient to filter the sample in a single operation, to avoid air intrusion and contamination.
- Avoid using glass syringes, as they may break due to filtration pressure. Depending on clogging, the applied pressure may exceed the maximum operating pressure; especially when using small syringes (1–2 mL), avoid applying excessive force.
- If filtration becomes difficult due to increased resistance, stop using the filter immediately.
- Always wear protective gloves and protective goggles when handling solvents and samples.

- If any contamination is detected, please perform a pre-rinse. The recommended pre-rinse volumes are as follows.

Filter Size	Pre-Rinse Volume
4 mm	0.5 - 1 mL
13 mm	3 - 5 mL
25 mm	10 - 20 mL

Sample considerations

For information regarding the effects of sample solutions on the filter, please refer to the Chemical Compatibility Chart. Due to variability in temperature, concentrations, duration of exposure, and other factors beyond our control, the information in the Chemical Compatibility Chart is provided without warranty, expressed or implied.

Disposal Precautions

After chemical solutions have been passed through this product, please dispose of it separately from general industrial waste and household waste, and in accordance with the laws and regulations of your local area.

Technical Support

If you encounter any performance defects, please contact your local Shimadzu sales office for assistance. However, please note that we cannot accept any claims when performance has degraded due to non-compliance with the operation procedures outlined above, or as a result of normal aging.

Chemical Compatibility Chart

Category	Sample solution	For IC
Acid	Acetic acid 10%	✓
	Acetic acid 30%	✓
	Hydrochloric acid	
	Nitric acid	
	Sulfuric acid	
Alkali	1.0 mol/L Potassium hydroxide (KOH)	✓
	3.0 mol/L Potassium hydroxide (KOH)	✓
	1.0 mol/L Sodium hydroxide (NaOH)	✓
	3.0 mol/L Sodium hydroxide (NaOH)	✓
Alcohol	Methanol	✓
	Ethanol	✓
	1-Propanol (n-propanol)	✓
	Isopropanol (2-propanol)	✓
	1-Butanol (n-butanol)	✓
	1-Pentanol (n-pentanol)	✓
	Ethylene glycol (ethane-1,2-diol)	✓
	Propylene glycol (1,2-propanediol)	✓
Glycerol (glycerin)	✓	
Others	Acetonitrile	
	N,N-Dimethylformamide (DMF)	
	Dimethyl sulfoxide (DMSO)	

This sheet and the appearance of the product may change without prior notice.

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