

# Multi-class pesticides analysis in challenging vegetable matrices using fast 5 msec MRM with 15 msec polarity switching

**IMSC 2012** PTh-176

Yuka Fujito<sup>1</sup>, Yusuke Inohana<sup>2</sup>, Kiyomi Arakawa<sup>2</sup>,  
Ichiro Hirano<sup>2</sup>

<sup>1</sup>Shimadzu Analytical & Measuring Center, Inc.,  
Kyoto, Japan ;

<sup>2</sup>Shimadzu Corporation, Kyoto, Japan

# Multi-class pesticides analysis in challenging vegetable matrices using fast 5 msec MRM with 15 msec polarity switching

## Introduction

Many regulatory authorities have established multi-class residual pesticides methods for the analysis of vegetables, fruits and other food stuffs. There is, however no global agreement on the provision of a target list of pesticides and this presents a risk with products moving between different regulatory requirements. In order to eliminate this risk, food safety laboratories need to ideally screen as many compounds as possible in a single run which may reach

maximum residual limits (MRL); typically 10 ppb in food matrices.

In this study, we report the application of ultra-fast 5 msec MRM with 15 msec polarity switching for the analysis of 138 pesticides in vegetable matrices (72 and 66 compounds measured by LC-QqQ and GC-QqQ in the European Union Reference Laboratory (EURL) method) .

## Materials and Methods

### Sample Preparation (QuEChERS EU method)



#### Food sample

Vegetables	Origin
Paprika (Sweet pepper)	New Zealand
Leek (Garlic chives)	Japan

#### Step 1 : Sample Extraction

1. Homogenize vegetables with food processor and homogenizer
2. Weigh 10 g homogenized sample
  - ← Add 10 mL acetonitrile
  - ← Add Salt-mixture\*1
3. Shake vigorously by hand 1 min.
4. Centrifuge for 5 min. at 4000 rpm (**Extract 1**)

- 4 g MgSO<sub>4</sub>
- 1 g NaCl
- 0.5 g Na<sub>2</sub>H citrate • 1.5H<sub>2</sub>O
- 1 g Na<sub>3</sub> citrate • 2H<sub>2</sub>O

\*1 : Citrate Extraction Tube (SIGMA ALDRICH)

#### Step 2 : Sample Clean up

5. Transfer 6 mL **Extract 1** into Dispersive SPE tube\*2
6. Shake vigorously by hand 2 min.
7. Centrifuge for 5 min. at 3000 rpm (**Extract 2**)
8. Transfer 1 mL **Extract 2** into a vial
9. Vortex sample to mix it
10. Filtrate sample with disposable filter

- containing
- 900 mg MgSO<sub>4</sub>
- 150 mg PSA
- 45 mg ENVI-Carb

\*2 : PSA/ENVI-Carb SPE Clean Up Tube 2 (SIGMA ALDRICH)

### LC/MS/MS analysis

#### Analytical Conditions

##### HPLC : Nexera UHPLC system

- Column : Shim-pack XR-ODSII (75 mm x 2 mmI.D., 2.2 μm)
- Mobile phase : A ; 2 mM ammonium formate containing 0.1% formic acid – water  
B ; Methanol
- Gradient program : 5% B (0-2.5 min.)→55% B (2.51-6 min.)→80% B (6.01-12 min.) →100% (12-15 min.)→5% (15.01-20 min.)
- Flow rate : 0.2 mL / min.
- Column temperature : 40°C

##### MS : LCMS-8040 Triple quadrupole mass spectrometer

- Ionization : ESI (Positive / Negative)
- Ion spray voltage : +4.5 kV / -3.5 kV
- MRM : 276 MRM transitions (2 MRMs / compound)  
Dwell time 5 msec. / Pause time 1 msec.

#### Features of LCMS-8040

- **5 times** higher sensitivity compared to LCMS-8030
- An ultra fast scan speed of **15000 u / sec.**
- An ultra fast polarity switching of **15 msec.**
- An ultra fast MRM transition speed of **555 ch./ sec.**



Fig. 1 LCMS-8040 Triple Quadrupole Mass Spectrometer

# Multi-class pesticides analysis in challenging vegetable matrices using fast 5 msec MRM with 15 msec polarity switching

## Results

### MRM of pesticide standards

Table 1 LOQs of 138 pesticides in the EURL method by LCMS-8040

Technique (on the EURL method)	LOQs < 10 ppb	LOQs > 10 ppb	Not Ionization
by LC-QqQ	<b>72 (100%)</b>	0 (0%)	0 (0%)
by GC-QqQ	<b>47 (71%)</b>	6 (9%)	13 (20%)

• 71 % of pesticides measured by GC-QqQ in the EURL method could achieved excellent LOQs (**0.08-10 ppb**).

#### Compounds for LC-QqQ

#### Compounds for GC-QqQ

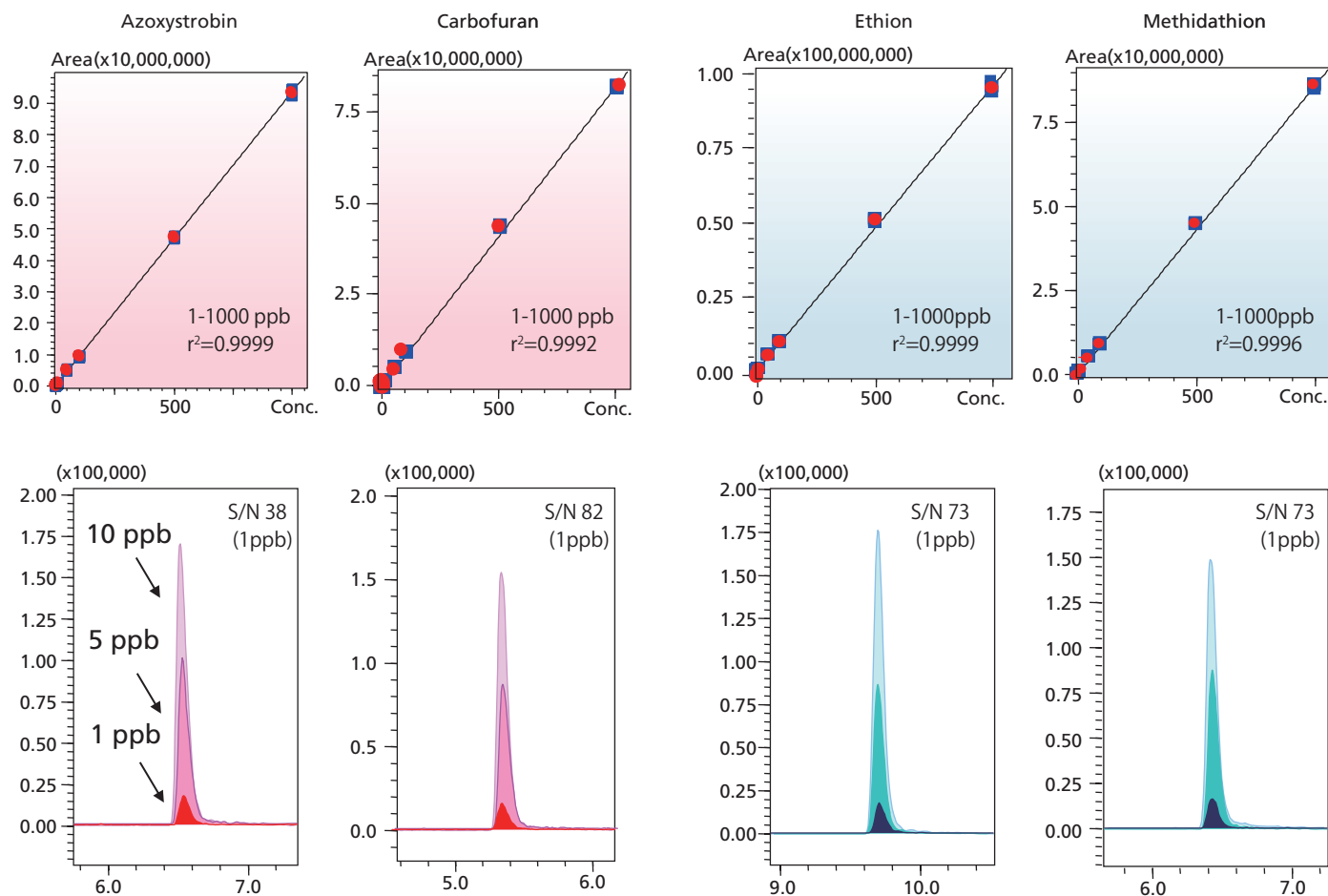


Fig. 2 Calibration curves and MRM chromatograms of typical pesticides

# Multi-class pesticides analysis in challenging vegetable matrices using fast 5 msec MRM with 15 msec polarity switching

## Recovery of pesticides in vegetable matrices


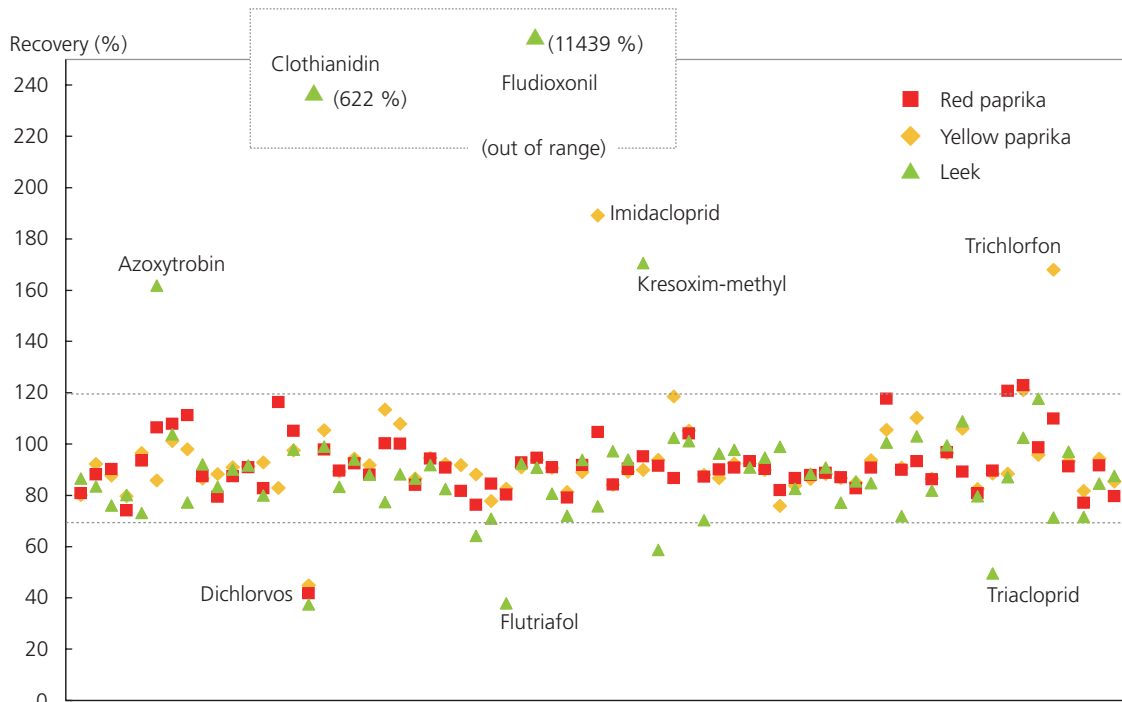


Fig. 3 Recovery of pesticides in the vegetable matrices (5 ppb spiked)

- Approximately **90%** of pesticides represented good recoveries in the range of 70-120% in all studied matrices.
- Some pesticides indicating up to 120% of recovery were also detected in the matrix blank (Fig. 4).

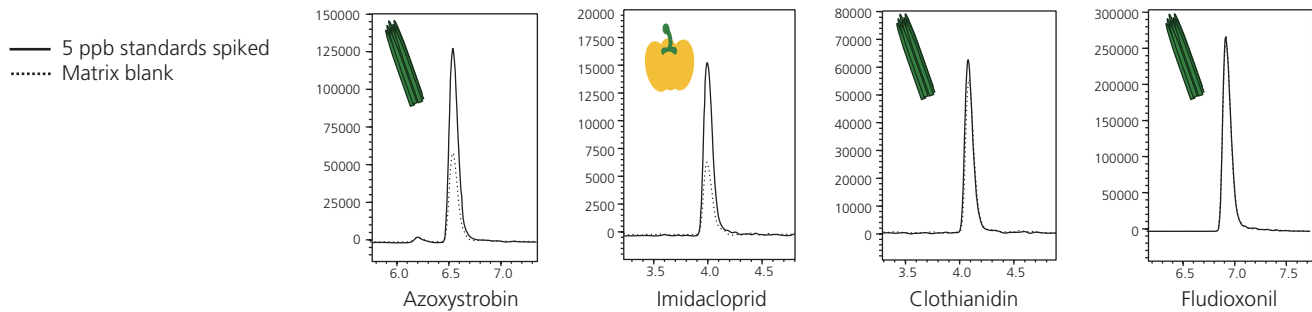

	Azoxystrobin	Imidacloprid	Clothianidin	Fludioxonil
5ppb standards spiked	8.19 ppb	8.99 ppb	25.78 ppb	556.58 ppb
Matrix blank	3.77 ppb	3.19 ppb	21.39 ppb	550.92 ppb
MRL (Japan)	5 ppm	3 ppm	15 ppm	10 ppm

Fig. 4 MRM Chromatograms and quantitative results of 4 pesticides

## Conclusion

- MRL of pesticide in vegetable matrices pre-treatment by QuEChERS method could be measured successfully using fast 5 msec MRM with 15 msec polarity switching.
- Majority of pesticides being suggested to GC-QqQ technology on the EURL method was successfully covered by LC-QqQ.



Shimadzu Corporation  
[www.shimadzu.com/an/](http://www.shimadzu.com/an/)

For Research Use Only. Not for use in diagnostic procedures.  
The content of this publication shall not be reproduced, altered or sold for any commercial purpose without the written approval of Shimadzu. The information contained herein is provided to you "as is" without warranty of any kind including without limitation warranties as to its accuracy or completeness. Shimadzu does not assume any responsibility or liability for any damage, whether direct or indirect, relating to the use of this publication. This publication is based upon the information available to Shimadzu on or before the date of publication, and subject to change without notice.